## CC100 Colony counter

*Please read the User Manual carefully before use, and follow all operating and safety instructions!*

## user manual

english, español

# User Manual

## LBX CC100 Colony Counter

### **Preface**

Users should read this Manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

### **Service**

If help is needed, you can always contact your dealer or Labbox via [www.labbox.com](http://www.labbox.com) (declare an incidence)

Please, provide the customer service representative with the following information:

- Serial number (on the back side)
- Description of the problem
- Your contact information

### **Warranty**

This instrument is guaranteed to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claim under the warranty, please contact your supplier.

# Contents

|   |                                      |
|---|--------------------------------------|
| Preface .....                                       | <b>¡Error! Marcador no definido.</b> |
| Service .....                                       | <b>¡Error! Marcador no definido.</b> |
| Warranty .....                                      | <b>¡Error! Marcador no definido.</b> |
| Contents .....                                      | 3                                    |
| 1.Safety Instructions.....                          | 4                                    |
| 2.Inspection .....                                  | 4                                    |
| 2.1 Receiving Inspection .....                      | 4                                    |
| 2.2 Listing of Items.....                           | 4                                    |
| 3.Assembly .....                                    | 5                                    |
| 4.Main part diagram .....                           | 6                                    |
| 5.Control panel .....                               | 7                                    |
| 6.Installation.....                                 | 7                                    |
| 7.Operation .....                                   | 8                                    |
| 8.Maintenance .....                                 | 10                                   |
| 9.Troubleshooting.....                              | 11                                   |
| 10.Software Instructions (Optional).....            | 12                                   |
| 11.Software Installation Guide.....                 | 12                                   |
| 12.Software operation guide – Data access .....     | 15                                   |
| 13.Software operation guide – Data deletion.....    | 18                                   |
| 14.Software operation guide – Data Query .....      | 20                                   |
| 15.Software operation guide – Data Print .....      | 21                                   |
| 16.Software operation guide – Data Saving .....     | 24                                   |
| 17.Software operation guide – Change Language ..... | 25                                   |

# 1. Safety Instructions

-This colony counter has been designed for laboratory usage only.

-Please read this manual carefully before installing or operating this equipment. The instrument shall not be modified in any way. Any modification will void the warranty and may result in potential safety hazard. We are not responsible for any injury or damage caused by modifying the instrument without authorization.

-To avoid the instrument from being burnt or disordered on operation, please use the power adaptor and USB connecting cable supplied by the manufacturer.

-Don't scratch the surface of the Petri dish holder to avoid affecting the illumination of the backlight.

-If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired. The instrument shall not be modified or altered in any way. Any modification or alteration will void the warranty, void the regulatory certifications and create a potential safety hazard.

-We are not responsible for any injury or damage caused by using the instrument for any non-intended purpose, or modifying the instrument by any person who is not authorized.

-Operating environment: Indoor use, 5 °C to 40 °C max. 80% RH max. Main supply voltage fluctuates up to  $\pm 10\%$  of the nominal voltage. Overvoltage category I. Pollution degree II.

## 2. Inspection

### 2.1 Receiving Inspection

Unpack the equipment carefully and check for any damages which may have arisen during transport. Please contact your supplier for technical support.

**Note:**

If there is any apparent damage to the system, please do not plug it into the power line.


### 2.2 Listing of Items

The package includes the following items:

| Items in CC100 Colony counter      | Qty |
|------------------------------------|-----|
| Colony counter                     | 1   |
| Magnifier                          | 1   |
| Magnifier holder                   | 1   |
| AC/DC switch adaptor               | 1   |
| Marking pen (black)                | 1   |
| Screwdriver                        | 1   |
| User Manual                        | 1   |
| Items in CC100 Software (Optional) | Qty |
| PC Software USB flash drive        | 1   |
| USB connecting cable               | 1   |
| Instruction manual                 | 1   |

Table 1


### 3. Assembly

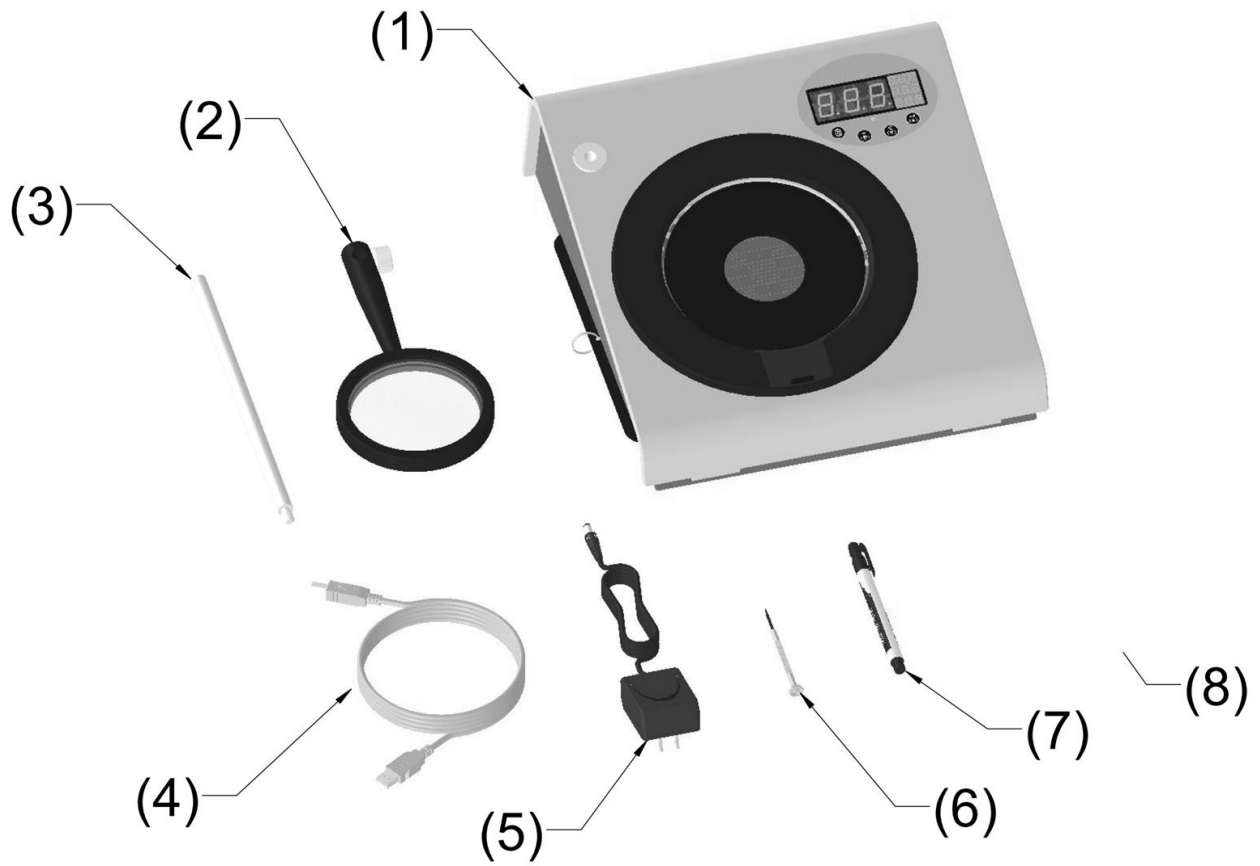| Item description                       | Number |
|--|--------|
| Colony counter                         | 1      |
| Magnifier                              | 2      |
| Magnifier holder                       | 3      |
| USB connecting cable (optional)        | 4      |
| AC/DC switching adaptor                | 5      |
| Screwdriver                            | 6      |
| Marking pen (black)                    | 7      |
| PC Software USB flash drive (optional) | 8      |

Table 2


#### 4. Main part diagram
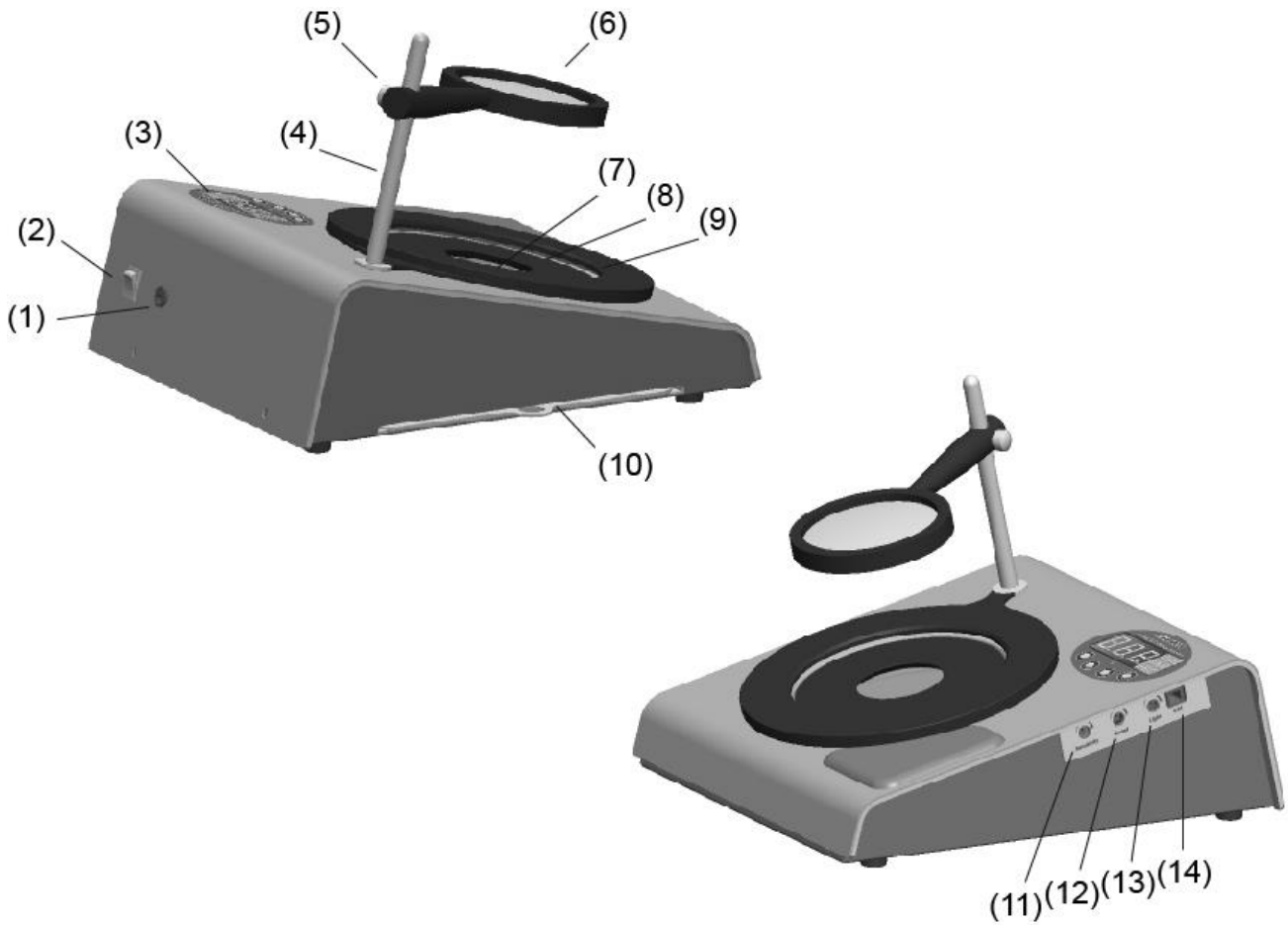
| Item description                         | Number |
|--|--------|
| Power socket                             | 1      |
| Power switch                             | 2      |
| Control panel                            | 3      |
| Magnifier holder                         | 4      |
| Adjustable screw                         | 5      |
| Magnifier                                | 6      |
| Petri dish position                      | 7      |
| Petri dish holder (60 mm)                | 8      |
| Adjustable petri dish holder (90~150 mm) | 9      |
| Background plate                         | 10     |
| Pressure sensitivity regulator           | 11     |
| Acoustic counting signal regulator       | 12     |
| Light intensity regulator                | 13     |
| USB connecting port                      | 14     |

Table 3


## 5. Control panel
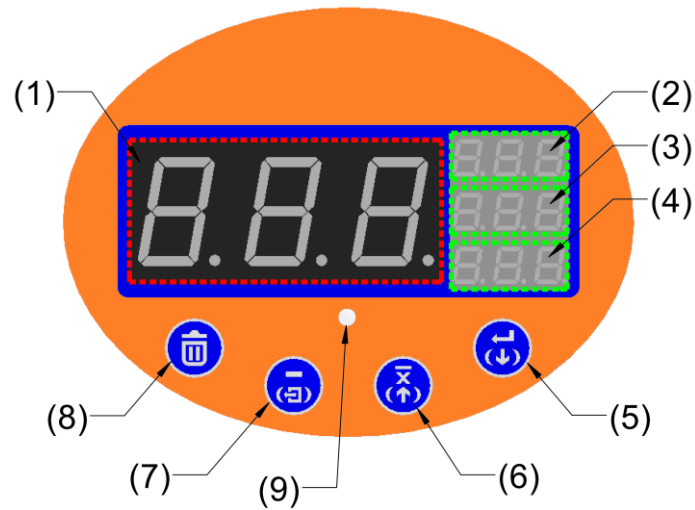
| Item description                                      | Number |
|---|--------|
| Counted value display (LED-A)                         | 1      |
| Storage value display (LED-B)                         | 2      |
| Storage value display (LED-C)                         | 3      |
| Storage value display (LED-D)                         | 4      |
| Save / Next   | 5      |
| Average / Previous                                    | 6      |
| Back / Exit   | 7      |
| Delete  | 8      |
| USB connecting indicator (green: connected, red: off) | 9      |

Table 4

## 6. Installation
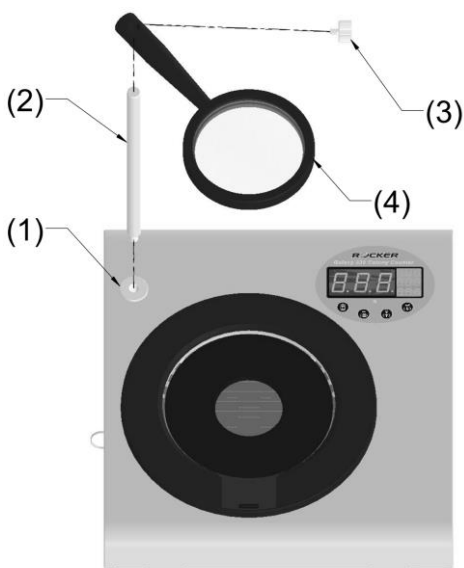
| Item description   | Number |
|--------------------|--------|
| Magnifier position | 1      |
| Magnifier holder   | 2      |
| Adjustable screw   | 3      |
| Magnifier          | 4      |

Table 5

1. Plug the power cord into the main unit and make sure the supplied voltage is within affordable range of power adaptor.
2. Install the magnifier on the main unit by following above diagram and adjust the height of magnifier properly.
3. If you want to transfer the counting data to a computer, please use USB cable provided to connect PC and colony counter and start the operation of software. For the first time use, please install the software by referring to the file of "PC Software Installation and Operation" on USB flash drive.
4. Please use the screwdriver provided to adjust the pressure sensitivity, acoustic volume, illumination required.
5. Change background plate (White/Black) if necessary
6. Please keep the Petri dish position and the magnifier clean to assure good light transmission.

## 7. Operation

### Key and Signal

#### (1) Normal Mode

|   |         |  |
|---|---------|--|
|    | SAVE    | Store the count in the memory                        |
|    | AVERAGE | Average saved counts<br>< display value on (LED-A) > |
|   | BACK    | Decrease the value on (LED-A)<br>(alarm with a beep) |
|  | DELETE  | Delete the count<br>< alarm with long beep >         |

#### (2) View Mode


|   |          |   |
|---|----------|---|
|  | EXIT     | 1). Press and hold this key for 1 second to enter<br>< View Mode ><br>when the display on (LED-A) is "000".<br>2). Press this key again to exit the <View Mode> |
|  | NEXT     | Check the next record   |
|  | PREVIOUS | Check the previous record   |
|  | DELETE   | Delete selected record on display (LED-B/C/D)   |

#### (3) Indicator Light

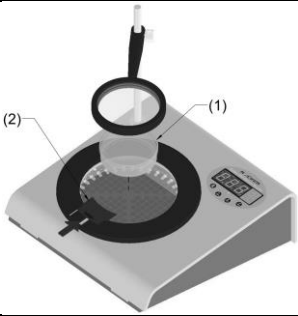- **Green light:** when the green light is ON means the connection is normal and the data transmission is successful. When light is flashing means the data is under transmission.
- **Red light:** when the red light is ON means the connection is abnormal and data transmission is failing.
- **No light:** it means the PC connection is not detectable.

## Operation under Normal Mode

|   |  |
|---|--|
|    | <p>Place the Petri dish (1) on the dish holder and slide the adjuster (2) properly.<br/>Attention: For the Petri dish with 60mm in diameter, you can use “Petri dish holder (60mm)” to fix the Petri dish.</p> |
|    | <p>Ensure the display on LED-A is "000" and no count display on LED-B、LED-C and LED-D before counting. If the count shows up, please press “DELETE” key to clear the memory.</p>                               |
|    | <p>Mark each colony with the marker pen. Every time a colony is marked, the apparatus will register the count on LED-A with a beep and counter advance.</p>  |
|   | <p>Press the “BACK” key if over counting occurs (alarm with a beep)</p>  |
|  | <p>After counting, press the “SAVE” key, the saved count will be displayed on LED-B.</p>   |
|  | <p>Place the next Petri dish and press &lt;SAVE &gt; key to resume the count. Repeat until all dishes have been counted. The saved counts will display on LED-B, LED-C, LED-D in order.</p>                    |
|  | <p>After you finish counting, press and hold “EXIT” key for 1 second to enter &lt; View Mode &gt; to check the saved counts. Press “EXIT” key again to be back to &lt;Normal Mode&gt;</p>                      |
|  | <p>Press the &lt;AVERAGE&gt; key to display the average count that calculated from all previously saved counts.</p>  |

## Notice:

- If the counted Petri dishes are over 3 batches, the display will only show the record of latest 3 batches. For the record of previous counts, it will be saved, but not shown on display.
- The main unit can store up to 100 records in memory. It will fail to work if the record is over 100. Please delete the record or activate "AVERAGE" facility, so the main unit can resume working.
- If the main unit connects to PC, the saved counts and average counts will be sent to the software on PC when you press "AVERAGE" key.
- Please refer to the file of "Installation and Operation of Software" on USB flash drive for operation of software.

### Operation under View Mode

|   |   |
|---|---|
|    | <p>Press and hold "EXIT" key for 1 second to enter &lt; View Mode &gt;.<br/>Notice: The display LED-A will show the total record numbers before you release the "EXIT" key.</p> |
|   | <p>After releasing the "EXIT" key, the saved counts will display on LED-B/C/D. The selected count will flash and LED-A will show the record number of selected count.</p>       |
|  | <p>Press the "NEXT" key to check the next record.<br/>It will alarm with a beep when it goes to the last record.</p>  |
|  | <p>Press the "PREVIOUS" key to check the previous record.<br/>It will alarm with a beep when it goes to the first record.</p>   |
|  | <p>Press "DELETE" key to delete flashed record. The alarm will sound with a long beep.</p>  |
|  | <p>Press "EXIT" key to be back to &lt;Normal Mode&gt;</p>   |

## 8. Maintenance

- After every use, please wipe the Petri dish position, the magnifier and surface of colony counter by 75% alcohol for sterilization and to gain good illumination.
- Please keep the instrument in well packed box to avoid dust if is not going to be used plan not to use it for a long time.

## 9. Troubleshooting

| Symptoms                                      | Possible causes and Solution  |
|---|---|
| <b>Device fails to start</b>                  | 1. Check the power supply conform to power specified (DC12V) on main unit<br>2. Check if the power switch is in the ON position<br>3. Malfunction of internal parts or internal tubing are loose → contact your distributor for repair      |
| <b>Poor LED lighting</b>                      | 1. Adjust the light intensity with a screwdriver<br>2. The backlight plate or mainboard is broken → contact your distributor for repair.  |
| <b>Acoustic counting signal is too low</b>    | 1. Adjust the volume with a screwdriver<br>2. Mainboard is broken → contact your distributor for repair.  |
| <b>Poor pressure sensitivity</b>              | 1. Adjust the sensitivity with a screwdriver<br>2. Mainboard is broken → contact your distributor for repair.   |
| <b>The main unit fails to connect with PC</b> | 1. Check if the USB cable is connected to the main unit and computer.<br>2. Check if the USB driver programming is installed<br>3. Make sure if the PC software is active.<br>4. Mainboard is broken → contact your distributor for repair. |

## 10. Software Instructions (Optional)

- This software can only be installed on PC or Laptop running Windows operating system. Please use the USB 2.0 cable supplied with the software package to ensure accurate data transfer.
- Recommended System Requirements:


|                  |   |
|------------------|---|
| Operating System | Windows XP SP3 32-bit / XP SP2 64-bit<br>Windows Vista SP2 32-bit / Vista SP2 64-bit<br>Windows 7 SP1 32-bit / 7 SP1 64-bit<br>Windows 8 32-bit / 8 64-bit ; Windows 8.1 32-bit / 8.1 64-bit<br>Windows 10 32-bit / 10 64-bit |
| CPU              | Pentium 4 1.5GHz above  |
| RAM              | 512MB above (1GB preferred)   |
| Hard Disk Space  | Installation 100MB above  |
| Interface        | USB Port  |
| Resolution       | 1024 × 768 Pixels   |

### Connection Status:


- When connected to the computer, the main screen displays  and the GREEN indicator on the main unit lights up.
- When disconnected of the computer, the main screen displays  and the RED indicator on the main unit lights up.

## 11. Software Installation Guide

- Insert the USB flash disk into one of the USB ports on the computer.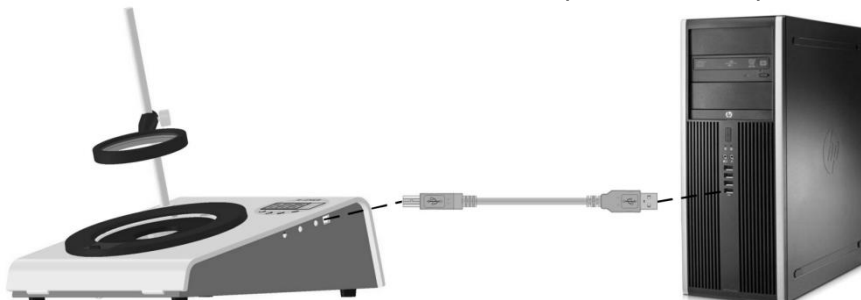

- Execute flash drive : \Galaxy-Setup\Galaxy330Install.exe
- Please click “Allow” or “Yes” if a prompt appears on screen asking for your permission.

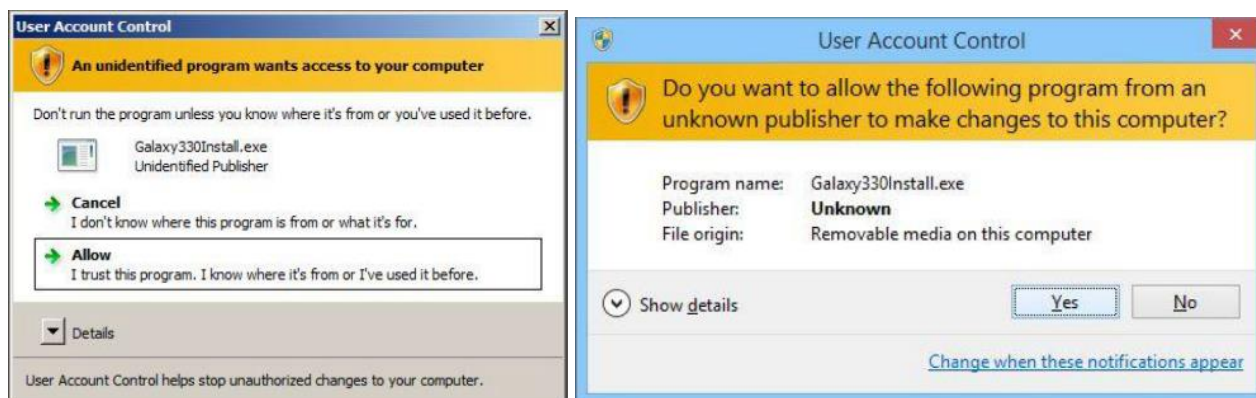
- Click “Setup” to install driver before software installation.  
(System auto detect 32- bit or 64 -bit version)


- Note: You must be logged in as an administrator to install this program.

**Driver Installation**

1. Click “Next” to start installation.
2. Click “I accept” and then click “Next” to begin installation.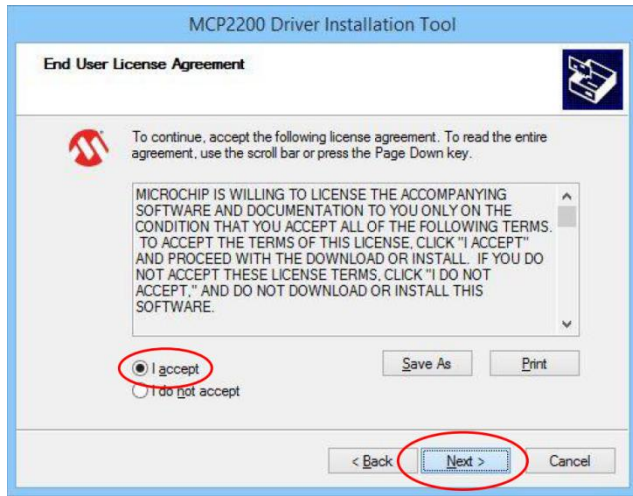

3. The driver installation begins and click “Install” to continue installation process.

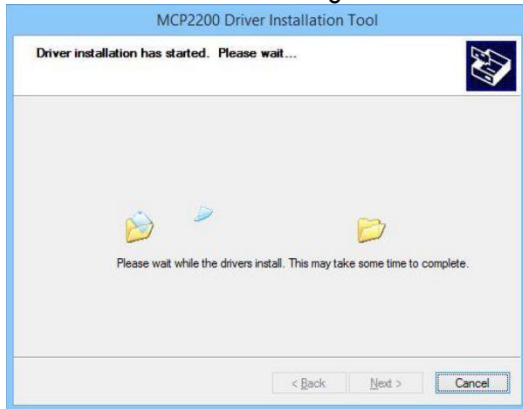4. Click "Finish" to complete the installation then begin installing software.


## Software Installation

1. Click "Next" to begin software installation.
2. Select installation folder then click "Next" to continue installation process.

3. Click "Install" to begin software installation.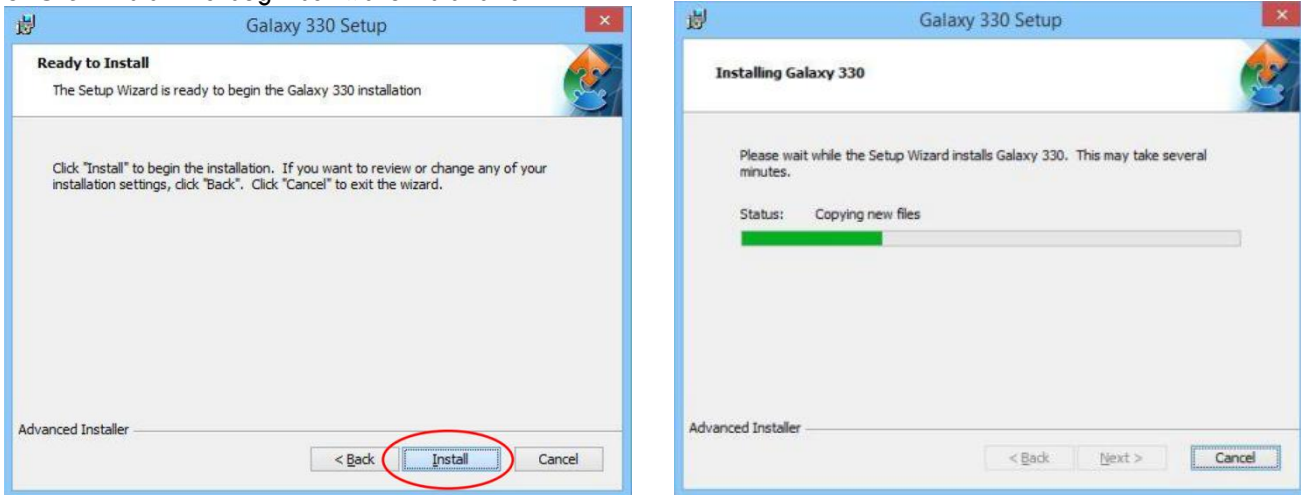

4. Click "Finish" to complete the installation

## 12. Software operation guide – Data access

### (1) Start Program

1. Click the Galaxy\_330 file icon (as shown below) under the Program Files folder or on the desktop.

2. Execute Galaxy 330 program then the main program screen will be displayed as the right figure.

**Note:** The main window display may vary depending on the operating system (OS), the user can set the main window properties by performing the following steps:




1. For Windows®, change window properties to Windows classic style.
2. For Windows® XP, click Control Panel → Display Properties → Appearance tab then change the color scheme to Windows classic.
3. For Windows® Vista / 7 / 8 / 8.1 / 10, click Control Panel → Personalization → Theme then change to the Windows Classic theme.

### (2) Input Sample data

Sample No: Input Sample Number  
Operator: Input Operator Name  
Customer: Input Customer Name  
Strain: Input Strain Name  
Media: Input Media Name  
Factor Scal.: Input Dilution Ratio  
Dilution Factor: Input Dilution Factor

### (3) Connect to Computer
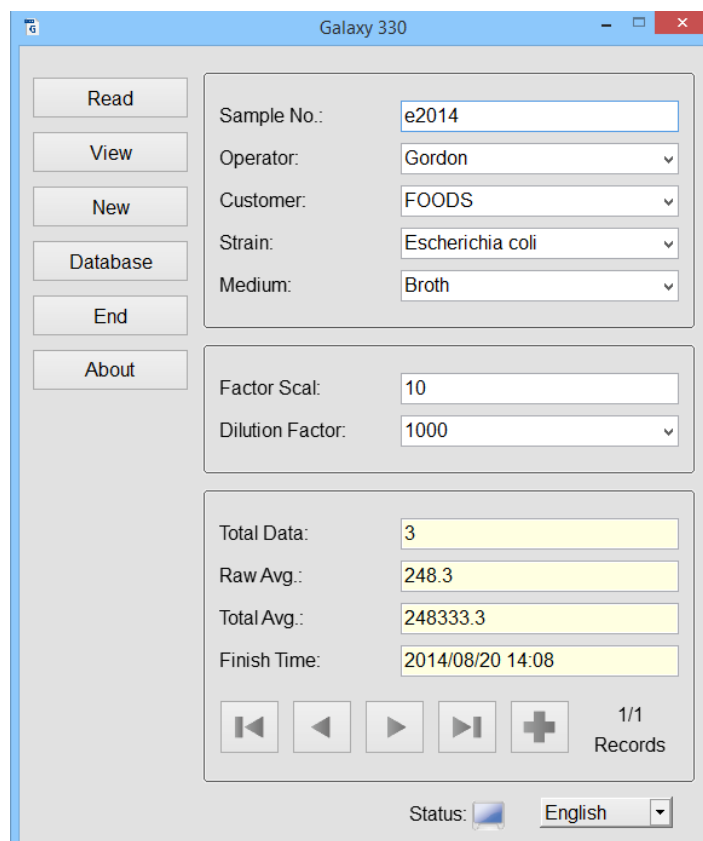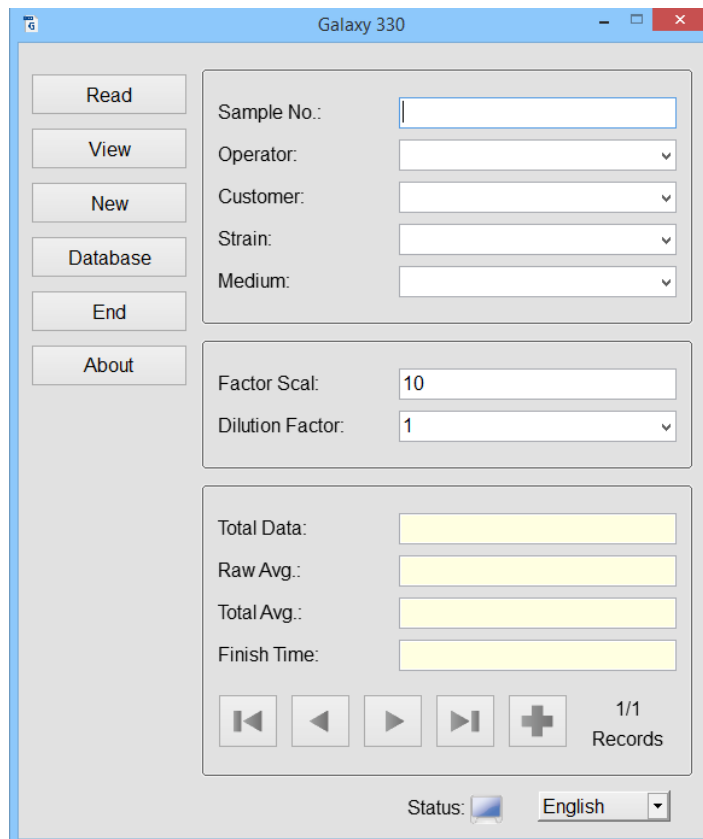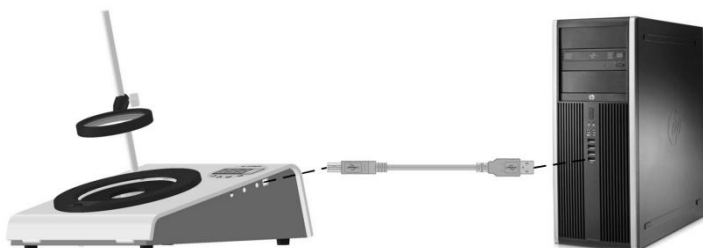
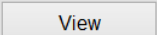
1. Connect to computer with USB cable

- Switch on Colony Counter then the Status indicator will display connection signal.
- When connected, the status indicator will display  and Colony Counter indicator will display "Green".
- When disconnected, the status indicator will display  and Colony Counter indicator will display "Red".

**Note: When the USB is connected, Windows® will automatically detect the hardware and perform driver, please be patient while in process.**

#### (4) Reading and Querying Data

- After counting bacterial colonies in the same sample and same dilution factor on each plate at Colony Counter, press "Average Button"  then each average count value will be automatically transmitted to the computer.
- Press  the main program screen and query all data received.

Galaxy 330 / View

|  |   |                                       |
|--|---|---------------------------------------|
| Sample No.: <input type="text" value="e2014"/><br>Operator: <input type="text" value="Gordon"/><br>Customer: <input type="text" value="FOODS"/><br>Strain: <input type="text" value="Escherichia coli"/> | Total Avg.: <input type="text" value="248333.3"/><br>Total Data: <input type="text" value="3"/> | <input type="button" value="Update"/> |
|--|---|---------------------------------------|

| Item | Operator | Strain           | Sample No. | Medium | Dilution Factor | Raw Count | Total Count            | Date                | Check |
|------|----------|------------------|------------|--------|-----------------|-----------|------------------------|---------------------|-------|
| 1    | Gordon   | Escherichia coli | e2014-001  | Broth  | 1000            | 243       | 2.43 × 10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 2    | Gordon   | Escherichia coli | e2014-002  | Broth  | 1000            | 253       | 2.53 × 10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 3    | Gordon   | Escherichia coli | e2014-003  | Broth  | 1000            | 249       | 2.49 × 10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |

- Press "EXIT" to return to the main program screen.

**Note: a set of data can contain max. 100 records and the number from 001 to 999 will be automatically added on counting data for different plate.**

**Note: in case the field cannot display all complete information, "... " will appear at the end of the field, then move mouse to the field, all complete information will be displayed after one second.**

## (5) Reading a new data

1. After completion of counting the following plates of bacterial colonies in the same sample and dilution factor at Colony Counter, press "Average Button"  then the next average counting data will be automatically transmitted to the computer again.
2. When turning into 2/2 Records, press each of the four keys below to switch display on counting data in different dilution factors up to 10 sets of data.

 To the first data

 To the previous data

 To the next data

 To the last data

Total Data:

Raw Avg.:

Total Avg.:

Finish Time:

2/2  
Records

3. Press  the main program screen to query all data received.

Galaxy 330 / View

Sample No.:

Operator:

Customer:

Strain:

Total Avg.:

Total Data:

| Item | Operator | Strain           | Sample No. | Medium | Dilution Factor | Raw Count | Total Count            | Date             | Check |
|------|----------|------------------|------------|--------|-----------------|-----------|------------------------|------------------|-------|
| 1    | Gordon   | Escherichia coli | e2014-001  | Broth  | 1000            | 243       | 2.43 × 10 <sup>5</sup> | 2014/08/20 14:08 | V     |
| 2    | Gordon   | Escherichia coli | e2014-002  | Broth  | 1000            | 253       | 2.53 × 10 <sup>5</sup> | 2014/08/20 14:08 | V     |
| 3    | Gordon   | Escherichia coli | e2014-003  | Broth  | 1000            | 249       | 2.49 × 10 <sup>5</sup> | 2014/08/20 14:08 | V     |
| 4    | Gordon   | Escherichia coli | e2014-004  | Broth  | 10000           | 24        | 2.40 × 10 <sup>5</sup> | 2014/08/20 14:15 | V     |
| 5    | Gordon   | Escherichia coli | e2014-005  | Broth  | 10000           | 25        | 2.50 × 10 <sup>5</sup> | 2014/08/20 14:15 | V     |
| 6    | Gordon   | Escherichia coli | e2014-006  | Broth  | 10000           | 29        | 2.90 × 10 <sup>5</sup> | 2014/08/20 14:15 | V     |

Press "EXIT" to return to the main program screen.

## (6) Viewing Data

Please repeat the previous steps for counting the same sample in different dilution factors. After completion of counting all different dilution factors for the same sample, press "View" to view all counting data, delete data, print data or save data as Excel file. Press "Exit" to return to the main program screen.

<<Please refer to respective instruction on how to delete, print or save data below>>

**Note:** in case the field cannot display all the information, "... " will appear at the end of the field, then move mouse to the field, all complete information will be displayed after one second.

## (7) Completing the count

After completing the counts of the previous sample, please press  and follow the above (1) ~ (6) steps to count dilution factors for the next sample.

**Note:**

a. After pressing the "New" key, a confirmation screen will appear. Press confirmation key to store all received data in database and then press  to view the set of data just saved;

b. This software can store up to max. 100,000 sets of data, once the number of the stored data exceeds the max. number, it will automatically delete the earliest data and then store the newly created data.

## 13. Software operation guide – Data deletion

### 1- Enter View Screen

During connection process, press  to query all received data in total 10 fields.

**Note: in case the field cannot display all complete information, "..." will appear at the end of the field, then move mouse to the field, all complete information will be displayed after one second.**

| Item | Operator | Strain           | Sample No. | Medium          | Dilution Factor | Raw Count | Total Count           | Date                | Check |
|------|----------|------------------|------------|-----------------|-----------------|-----------|-----------------------|---------------------|-------|
| 1    | Gordon   | Escherichia coli | e2014-001  | Synthetic me... | 1000            | 243       | 2.43 ×10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 2    | Gordon   | Escherichia coli | e2014-002  | Synthetic me... | 1000            | 253       | 2.53 ×10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 3    | Gordon   | Escherichia coli | e2014-003  | Synthetic me... | 1000            | 249       | 2.49 ×10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 4    | Gordon   | Escherichia coli | e2014-004  | Synthetic me... | 10000           | 24        | 2.40 ×10 <sup>5</sup> | 2014/08/20<br>14:15 | V     |
| 5    | Gordon   | Escherichia coli | e2014-005  | Synthetic me... | 10000           | 25        | 2.50 ×10 <sup>5</sup> | 2014/08/20<br>14:15 | V     |
| 6    | Gordon   | Escherichia coli | e2014-006  | Synthetic me... | 10000           | 29        | 2.90 ×10 <sup>5</sup> | 2014/08/20<br>14:15 | V     |

Synthetic media

|                 |   |
|-----------------|---|
| Item            | Data Number                               |
| Operator        | Operator Name                             |
| Strain          | Strain Name                               |
| Sample No       | Counted Sample Number                     |
| Media           | Media Name                                |
| Dilution Factor | Dilution Factor for a concentrated sample |
| Raw Count       | Colony count number                       |
| Total Count     | The number of colonies after conversion   |
| Date            | The data-receiving date                   |
| Check           | Check if the data is valid                |

### 2- Select the data you want to delete

If any incorrect count data is found, move mouse cursor to the Check field.

| Item | Operator | Strain           | Sample No. | Medium          | Dilution Factor | Raw Count | Total Count           | Date                | Check |
|------|----------|------------------|------------|-----------------|-----------------|-----------|-----------------------|---------------------|-------|
| 1    | Gordon   | Escherichia coli | e2014-001  | Synthetic me... | 1000            | 243       | 2.43 ×10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 2    | Gordon   | Escherichia coli | e2014-002  | Synthetic me... | 1000            | 253       | 2.53 ×10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 3    | Gordon   | Escherichia coli | e2014-003  | Synthetic me... | 1000            | 249       | 2.49 ×10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 4    | Gordon   | Escherichia coli | e2014-004  | Synthetic me... | 10000           | 24        | 2.40 ×10 <sup>5</sup> | 2014/08/20<br>14:15 | V     |
| 5    | Gordon   | Escherichia coli | e2014-005  | Synthetic me... | 10000           | 25        | 2.50 ×10 <sup>5</sup> | 2014/08/20<br>14:15 | V     |
| 6    | Gordon   | Escherichia coli | e2014-006  | Synthetic me... | 10000           | 29        | 2.90 ×10 <sup>5</sup> | 2014/08/20<br>14:15 | V     |

Synthetic media

Click the left mouse button to "V" sign to make the field become blank

| Item | Operator | Strain           | Sample No. | Medium          | Dilution Factor | Raw Count | Total Count           | Date                | Check |
|------|----------|------------------|------------|-----------------|-----------------|-----------|-----------------------|---------------------|-------|
| 1    | Gordon   | Escherichia coli | e2014-001  | Synthetic me... | 1000            | 243       | 2.43 ×10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 2    | Gordon   | Escherichia coli | e2014-002  | Synthetic me... | 1000            | 253       | 2.53 ×10 <sup>5</sup> | 2014/08/20<br>14:08 |       |
| 3    | Gordon   | Escherichia coli | e2014-003  | Synthetic me... | 1000            | 249       | 2.49 ×10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 4    | Gordon   | Escherichia coli | e2014-004  | Synthetic me... | 10000           | 24        | 2.40 ×10 <sup>5</sup> | 2014/08/20<br>14:15 | V     |
| 5    | Gordon   | Escherichia coli | e2014-005  | Synthetic me... | 10000           | 25        | 2.50 ×10 <sup>5</sup> | 2014/08/20<br>14:15 | V     |
| 6    | Gordon   | Escherichia coli | e2014-006  | Synthetic me... | 10000           | 29        | 2.90 ×10 <sup>5</sup> | 2014/08/20<br>14:15 |       |

**Note: The blank field indicates the data will be deleted and re-click on the blank field will show the following screen. Click on the unwanted data and press "Update" to confirm the deletion.**

|             |   |             |                                       |                                       |
|-------------|---|-------------|---------------------------------------|---------------------------------------|
| Sample No.: | <input type="text" value="e2014"/>            | Total Avg.: | <input type="text" value="245500.0"/> | <input type="button" value="Update"/> |
| Operator:   | <input type="text" value="Gordon"/>           | Total Data: | <input type="text" value="4"/>        |                                       |
| Customer:   | <input type="text" value="FOODS"/>            |             |                                       |                                       |
| Strain:     | <input type="text" value="Escherichia coli"/> |             |                                       |                                       |

Press "Yes" to reconfirm the data deletion. (unable to be recovered after deletion)

Galaxy 330 / Warning

Change Data!! Yes/No

Shown as below after data deletion.

Galaxy 330 / View

|             |   |             |                                       |                                       |
|-------------|---|-------------|---------------------------------------|---------------------------------------|
| Sample No.: | <input type="text" value="e2014"/>            | Total Avg.: | <input type="text" value="245500.0"/> | <input type="button" value="Update"/> |
| Operator:   | <input type="text" value="Gordon"/>           | Total Data: | <input type="text" value="4"/>        |                                       |
| Customer:   | <input type="text" value="FOODS"/>            |             |                                       |                                       |
| Strain:     | <input type="text" value="Escherichia coli"/> |             |                                       |                                       |

| Item | Operator | Strain           | Sample No. | Medium          | Dilution Factor | Raw Count | Total Count           | Date                | Check |
|------|----------|------------------|------------|-----------------|-----------------|-----------|-----------------------|---------------------|-------|
| 1    | Gordon   | Escherichia coli | e2014-001  | Synthetic me... | 1000            | 243       | 2.43 ×10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 2    | Gordon   | Escherichia coli | e2014-003  | Synthetic me... | 1000            | 249       | 2.49 ×10 <sup>5</sup> | 2014/08/20<br>14:08 | V     |
| 3    | Gordon   | Escherichia coli | e2014-004  | Synthetic me... | 10000           | 24        | 2.40 ×10 <sup>5</sup> | 2014/08/20<br>14:15 | V     |
| 4    | Gordon   | Escherichia coli | e2014-005  | Synthetic me... | 10000           | 25        | 2.50 ×10 <sup>5</sup> | 2014/08/20<br>14:15 | V     |

(3) After completion of data deletion and press "Exit" to return to the main program screen.

# 14. Software operation guide – Data Query

-Enter View Screen

Press "Database" in the main program screen to enter view screen.

The screenshot shows the 'Galaxy 330 / Database' window. It features a search criteria section with the following fields:

- Sample No.:
- Operator:
- Customer:
- Strain:
- Start Date:
- End Date:
- Total Data:
- Total Avg.:

Buttons for 'Search' and 'Clear' are located to the right of the date and total data fields. Below the search criteria is a table with the following columns: Item, Operator, Customer, Strain, Sample No., Medium, Dilution Factor, Raw Count, Total Count, and Date. To the right of the table are buttons for 'Title', 'Print', 'Save as', and 'Exit'.

The six fields below can be set as query conditions.

|            |
|------------|
| Sample No. |
| Operator   |
| Customer   |
| Strain     |
| Start Date |
| End Date   |

Note: The query condition can accept English capital, small letters or incomplete characters.

-Start Search

Press "Search" after inputting the query condition then all queried data will appear.

Press "Clear" to clear all displayed data before re-query.

The screenshot shows the 'Galaxy 330 / Database' window after a search. The search criteria are the same as in the previous screenshot. The 'Search' and 'Clear' buttons are highlighted with a red box. The table below shows the results of the search:

| Item | Operator | Customer | Strain          | Sample No. | Medium          | Dilution Factor | Raw Count | Total Count            | Date             |
|------|----------|----------|-----------------|------------|-----------------|-----------------|-----------|------------------------|------------------|
| 1    | Gordon   | FOODS    | Escherichia ... | e2014-001  | Synthetic me... | 1000            | 243       | 2.43 × 10 <sup>5</sup> | 2014/08/20 14:08 |
| 2    | Gordon   | FOODS    | Escherichia ... | e2014-003  | Synthetic me... | 1000            | 249       | 2.49 × 10 <sup>5</sup> | 2014/08/20 14:08 |
| 3    | Gordon   | FOODS    | Escherichia ... | e2014-004  | Synthetic me... | 10000           | 24        | 2.40 × 10 <sup>5</sup> | 2014/08/20 14:15 |
| 4    | Gordon   | FOODS    | Escherichia ... | e2014-005  | Synthetic me... | 10000           | 25        | 2.50 × 10 <sup>5</sup> | 2014/08/20 14:15 |

Note: This software can display up to 1000 sets of data. Please narrow your search to avoid excessive data.

## -Search Result

The search result will be displayed as above in total 10 fields incl. Item No., Operator, Customer, Strain, Sample No, Media, Dilution Factor, Raw Count, Total Count and Date.

**Note: in case the field cannot display all complete information, "..."** will appear at the end of the field, then move mouse to the field, all complete information will be displayed after one second.

## 15. Software operation guide – Data Print

### (1) Search the data for printing

Search the data for printing by pressing “View” or “Database” in the main program screen and a sub-screen will appear. Press “Print” and set the printer in portrait, A4 format then the data list will be printed out.

The screenshot shows the 'Galaxy 330 / View' software interface. At the top, there are input fields for 'Sample No.' (e2014), 'Operator' (Gordon), 'Customer' (FOODS), and 'Strain' (Escherichia coli). To the right, there are fields for 'Total Avg.: 245500.0' and 'Total Data: 4', with an 'Update' button. Below these is a table with 10 columns: Item, Operator, Strain, Sample No., Medium, Dilution Factor, Raw Count, Total Count, Date, and Check. The table contains 4 rows of data. To the right of the table is a vertical stack of buttons: 'Title', 'Print' (highlighted with a red box), 'Save as', and 'Exit'.

| Item | Operator | Strain           | Sample No. | Medium          | Dilution Factor | Raw Count | Total Count            | Date             | Check |
|------|----------|------------------|------------|-----------------|-----------------|-----------|------------------------|------------------|-------|
| 1    | Gordon   | Escherichia coli | e2014-001  | Synthetic me... | 1000            | 243       | 2.43 × 10 <sup>5</sup> | 2014/08/20 14:08 | V     |
| 2    | Gordon   | Escherichia coli | e2014-003  | Synthetic me... | 1000            | 249       | 2.49 × 10 <sup>5</sup> | 2014/08/20 14:08 | V     |
| 3    | Gordon   | Escherichia coli | e2014-004  | Synthetic me... | 10000           | 24        | 2.40 × 10 <sup>5</sup> | 2014/08/20 14:15 | V     |
| 4    | Gordon   | Escherichia coli | e2014-005  | Synthetic me... | 10000           | 25        | 2.50 × 10 <sup>5</sup> | 2014/08/20 14:15 | V     |

The screenshot shows the 'Print' dialog box. It has a 'Printer' section with 'Name: Microsoft XPS Document Writer' and a 'Properties...' button. Below that, it shows 'Status: Ready', 'Type: Microsoft XPS Document Writer v4', and 'Where: PORTPROMPT:'. There are also 'Print range' options: 'All' (selected), 'Pages' (with 'from:' and 'to:' fields), and 'Selection'. The 'Copies' section has 'Number of copies: 1' and a 'Collate' checkbox. At the bottom are 'OK' and 'Cancel' buttons.

# COLONY COUNTING REPORT

| Item          | Operator | Strain           | Sample No. | Medium          | Dilution Factor                 | Raw Count | Total Count           | Date             |
|---------------|----------|------------------|------------|-----------------|---------------------------------|-----------|-----------------------|------------------|
| 1             | Gordon   | Escherichia coli | e2014-001  | Synthetic me... | 1000                            | 243       | 2.43 ×10 <sup>5</sup> | 2014/08/20 14:08 |
| 2             | Gordon   | Escherichia coli | e2014-003  | Synthetic me... | 1000                            | 249       | 2.49 ×10 <sup>5</sup> | 2014/08/20 14:08 |
| 3             | Gordon   | Escherichia coli | e2014-004  | Synthetic me... | 10000                           | 24        | 2.40 ×10 <sup>5</sup> | 2014/08/20 14:15 |
| 4             | Gordon   | Escherichia coli | e2014-005  | Synthetic me... | 10000                           | 25        | 2.50 ×10 <sup>5</sup> | 2014/08/20 14:15 |
|               |          |                  |            |                 |                                 |           |                       |                  |
|               |          |                  |            |                 |                                 |           |                       |                  |
|               |          |                  |            |                 |                                 |           |                       |                  |
|               |          |                  |            |                 |                                 |           |                       |                  |
|               |          |                  |            |                 |                                 |           |                       |                  |
|               |          |                  |            |                 |                                 |           |                       |                  |
|               |          |                  |            |                 |                                 |           |                       |                  |
|               |          |                  |            |                 |                                 |           |                       |                  |
|               |          |                  |            |                 |                                 |           |                       |                  |
|               |          |                  |            |                 |                                 |           |                       |                  |
| Total data: 4 |          |                  |            |                 | Average: 2.455 ×10 <sup>5</sup> |           |                       |                  |

Print date: 2014/08/20 15:46

1

*Note: The printing format is designed for the portrait, A4 paper only so please don't use another size. To print the other paper formats. Please save the data as Excel file and edit the file with Excel before printing.*

**(2) Printing the report with title**

Press "Title" in the sub-screen from pressing "View" or "Database" to print the report with title.

Galaxy 330 / View

|   |   |                                       |
|---|---|---------------------------------------|
| Sample No.: <input type="text" value="e2014"/>        | Total Avg.: <input type="text" value="245500.0"/> | <input type="button" value="Update"/> |
| Operator: <input type="text" value="Gordon"/>         | Total Data: <input type="text" value="4"/>        |                                       |
| Customer: <input type="text" value="FOODS"/>          |   |                                       |
| Strain: <input type="text" value="Escherichia coli"/> |   |                                       |

| Item | Operator | Strain           | Sample No. | Medium          | Dilution Factor | Raw Count | Total Count           | Date             | Check |
|------|----------|------------------|------------|-----------------|-----------------|-----------|-----------------------|------------------|-------|
| 1    | Gordon   | Escherichia coli | e2014-001  | Synthetic me... | 1000            | 243       | 2.43 ×10 <sup>5</sup> | 2014/08/20 14:08 | V     |
| 2    | Gordon   | Escherichia coli | e2014-003  | Synthetic me... | 1000            | 249       | 2.49 ×10 <sup>5</sup> | 2014/08/20 14:08 | V     |
| 3    | Gordon   | Escherichia coli | e2014-004  | Synthetic me... | 10000           | 24        | 2.40 ×10 <sup>5</sup> | 2014/08/20 14:15 | V     |
| 4    | Gordon   | Escherichia coli | e2014-005  | Synthetic me... | 10000           | 25        | 2.50 ×10 <sup>5</sup> | 2014/08/20 14:15 | V     |

**ROCKER**

Fill in all required information after the following form appears.

Galaxy 330 / Title

Company:

Department: Sales Department

Tel:

Fax:

Address:

E-mail:

Operator: Gordon

Approval: Tim

Print Title

Submit

Cancel

Load


Save

Make sure that "Print Title" has been ticked after filling in all the title information.


Note: if "Print Title" is not ticked then the title won't be printed out.

#### (4) Saving Data

Press "Save" to save all title information.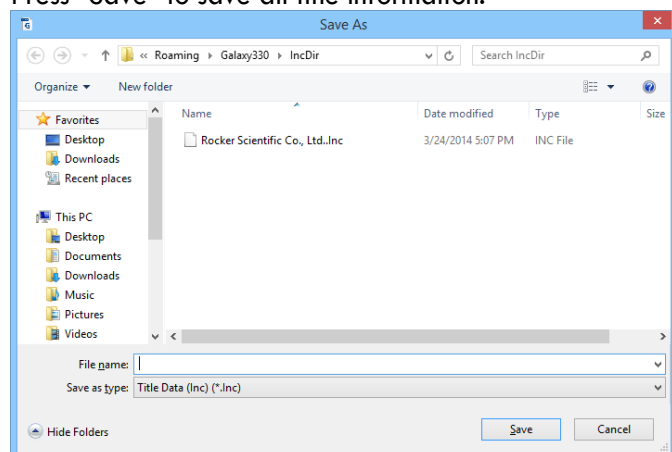

All saved title information can then be re-loaded back into the program by pressing "Load".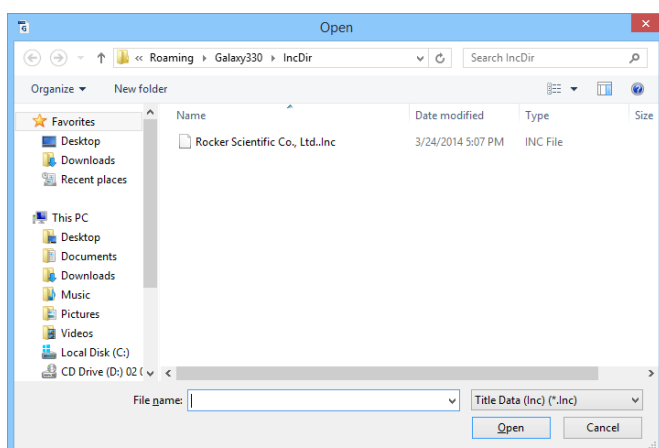

#### (5) Print the report with title

Ensure all title information are correct and press "Submit" to return to the previous screen. After pressing "Print" and selecting the printer, the report with title can be printed out.

#### Note:

- Press "Cancel" to abort any amended title information and all previous title information will keep unchanged.
- Uncheck "Print Title" if the title information is not required for the report.

COLONY COUNTING REPORT

Company:

Department: Sales Department

Address:

Tel:

Fax:

E-mail:

Operator: Gordon

Approval: Tim



| Item          | Operator | Strain           | Sample No. | Medium          | Dilution Factor                  | Raw Count | Total Count            | Date             |
|---------------|----------|------------------|------------|-----------------|----------------------------------|-----------|------------------------|------------------|
| 1             | Gordon   | Escherichia coli | e2014-001  | Synthetic me... | 1000                             | 243       | 2.43 × 10 <sup>5</sup> | 2014/08/20 14:08 |
| 2             | Gordon   | Escherichia coli | e2014-003  | Synthetic me... | 1000                             | 249       | 2.49 × 10 <sup>5</sup> | 2014/08/20 14:08 |
| 3             | Gordon   | Escherichia coli | e2014-004  | Synthetic me... | 10000                            | 24        | 2.40 × 10 <sup>5</sup> | 2014/08/20 14:15 |
| 4             | Gordon   | Escherichia coli | e2014-005  | Synthetic me... | 10000                            | 25        | 2.50 × 10 <sup>5</sup> | 2014/08/20 14:15 |
| Total data: 4 |          |                  |            |                 | Average: 2.455 × 10 <sup>5</sup> |           |                        |                  |

Print date: 2014/08/20 15:48

## 16. Software operation guide – Data Saving

(1) Search the data to be saved

Search the data to be saved through the sub-screen from pressing "View" or "Database".  
Click on "Save as" to save the data as CSV file of the Excel.

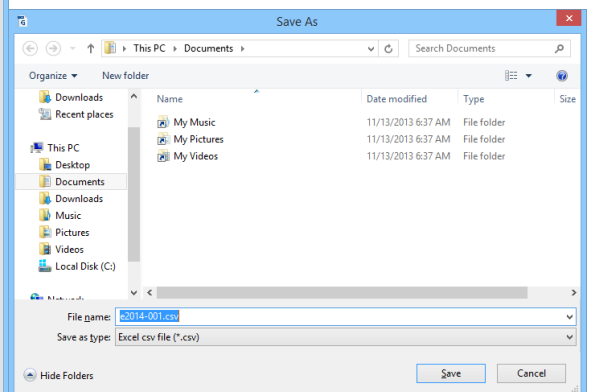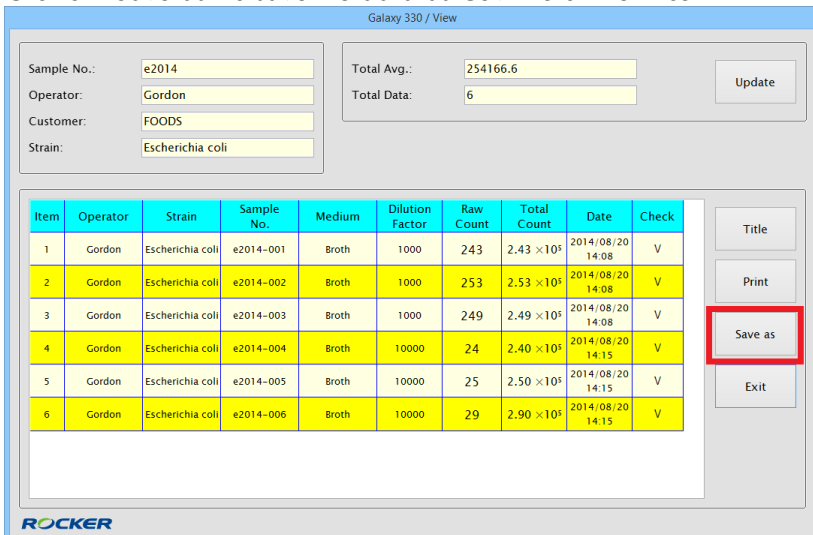(2) Open .CSV File with Excel

After opening Excel, select open the old file (.csv file).

| Item | Operator | Customer | Strain           | Sample No. | Medium          | Dilution Factor | Raw Count | Total Count | Date            |
|------|----------|----------|------------------|------------|-----------------|-----------------|-----------|-------------|-----------------|
| 12   | 1 Gordon | FOODS    | Escherichia coli | e2014-001  | Synthetic media | 1000            | 243       | 2.43E+05    | 2014/8/20 14:08 |
| 13   | 2 Gordon | FOODS    | Escherichia coli | e2014-003  | Synthetic media | 1000            | 249       | 2.49E+05    | 2014/8/20 14:08 |
| 14   | 3 Gordon | FOODS    | Escherichia coli | e2014-004  | Synthetic media | 10000           | 24        | 2.40E+05    | 2014/8/20 14:15 |
| 15   | 4 Gordon | FOODS    | Escherichia coli | e2014-005  | Synthetic media | 10000           | 25        | 2.50E+05    | 2014/8/20 14:15 |

After opening the .csv file, all the data can be tidied as desired.

**Note:**

- a. Whether ticking “Print Title” or not, all existing title information will be included in the .CSV file. However, these title information can be edited or deleted as requested.
- b. The data of total count will be expressed in scientific notation.

## 17. Software operation guide – Change Language

**CHANGE LANGUAGE:**

Move mouse cursor to the right button at the main program screen to switch English / Traditional Chinese/Simple Chinese lanaguage interfaces:

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The screenshot shows the Galaxy 330 / Database software interface. At the top, there are search criteria fields: Sample No., Operator, Customer, and Strain (all dropdown menus); Start Date and End Date (date pickers); Total Data (text input); and Total Avg. (text input). To the right of these fields are Search and Clear buttons. Below the search criteria is a large table with columns: Item, Operator, Customer, Strain, Sample No., Medium, Dilution Factor, Raw Count, Total Count, and Date. The table is currently empty. To the right of the table are buttons for Title, Print, Save as, and Exit.

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The screenshot shows the Galaxy 330 / Database software interface with search results. The search criteria fields are: Sample No. (empty), Operator (Gordon), Customer (empty), Strain (empty), Start Date (2014/08/20), End Date (2014/08/20), Total Data (4), and Total Avg. (245500.0). The Search and Clear buttons are highlighted with a red box. The data table now contains four rows of results:

| Item | Operator | Customer | Strain          | Sample No. | Medium          | Dilution Factor | Raw Count | Total Count            | Date             |
|------|----------|----------|-----------------|------------|-----------------|-----------------|-----------|------------------------|------------------|
| 1    | Gordon   | FOODS    | Escherichia ... | e2014-001  | Synthetic me... | 1000            | 243       | 2.43 × 10 <sup>5</sup> | 2014/08/20 14:08 |
| 2    | Gordon   | FOODS    | Escherichia ... | e2014-003  | Synthetic me... | 1000            | 249       | 2.49 × 10 <sup>5</sup> | 2014/08/20 14:08 |
| 3    | Gordon   | FOODS    | Escherichia ... | e2014-004  | Synthetic me... | 10000           | 24        | 2.40 × 10 <sup>5</sup> | 2014/08/20 14:15 |
| 4    | Gordon   | FOODS    | Escherichia ... | e2014-005  | Synthetic me... | 10000           | 25        | 2.50 × 10 <sup>5</sup> | 2014/08/20 14:15 |

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The screenshot shows the 'Galaxy 330 / View' software interface. It features a search form on the left with fields for Sample No., Operator, Customer, and Strain. On the right, there are summary statistics for Total Avg. and Total Data, along with an 'Update' button. Below the search form is a table with 10 columns: Item, Operator, Strain, Sample No., Medium, Dilution Factor, Raw Count, Total Count, Date, and Check. The table contains 4 rows of data. To the right of the table is a vertical menu with buttons for 'Title', 'Print', 'Save as', and 'Exit'. The 'Print' button is highlighted with a red rectangle.

| Item | Operator | Strain           | Sample No. | Medium          | Dilution Factor | Raw Count | Total Count            | Date             | Check |
|------|----------|------------------|------------|-----------------|-----------------|-----------|------------------------|------------------|-------|
| 1    | Gordon   | Escherichia coli | e2014-001  | Synthetic me... | 1000            | 243       | 2.43 × 10 <sup>5</sup> | 2014/08/20 14:08 | V     |
| 2    | Gordon   | Escherichia coli | e2014-003  | Synthetic me... | 1000            | 249       | 2.49 × 10 <sup>5</sup> | 2014/08/20 14:08 | V     |
| 3    | Gordon   | Escherichia coli | e2014-004  | Synthetic me... | 10000           | 24        | 2.40 × 10 <sup>5</sup> | 2014/08/20 14:15 | V     |
| 4    | Gordon   | Escherichia coli | e2014-005  | Synthetic me... | 10000           | 25        | 2.50 × 10 <sup>5</sup> | 2014/08/20 14:15 | V     |

The screenshot shows the 'Print' dialog box. It has a 'Printer' section with a dropdown menu set to 'Microsoft XPS Document Writer' and a 'Properties...' button. Below this, it shows the printer's status as 'Ready', type as 'Microsoft XPS Document Writer v4', and location as 'PORTPROMPT:'. There are also 'Print range' and 'Copies' sections. The 'Print range' section has radio buttons for 'All', 'Pages' (with 'from' and 'to' fields), and 'Selection'. The 'Copies' section has a 'Number of copies' spinner set to 1 and a 'Collate' checkbox. At the bottom are 'OK' and 'Cancel' buttons.

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