CC100 Colony counter

Please read the User Manual carefully before use, and follow all operating and safety instructions!

user manual

english, español

User Manual

LBX CC100 Colony Counter

Preface

Users should read this Manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

Service

If help is needed, you can always contact your dealer or Labbox via www.labbox.com (declare an incidence)

Please, provide the customer service representative with the following information:

- Serial number (on the back side)
- Description of the problem
- Your contact information

Warranty

This instrument is guaranteed to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claim under the warranty, please contact your supplier.

Contents

Preface	¡Error! Marcador no definido.
Service	¡Error! Marcador no definido.
Warranty	¡Error! Marcador no definido.
Contents	3
1.Safety Instructions.....	3
2.Inspection	4
2.1 Receiving Inspection	4
2.2 Listing of Items.....	4
3.Assembly	5
4.Main part diagram	6
5.Control panel	7
6.Installation.....	7
7.Operation	8
8.Maintenance	10
9.Troubleshooting.....	11
10.Software Instructions (Optional).....	12
11.Software Installation Guide.....	12
12.Software operation guide – Data access	15
13.Software operation guide – Data deletion.....	18
14.Software operation guide – Data Query	20
15.Software operation guide – Data Print	21
16.Software operation guide – Data Saving	24
17.Software operation guide – Change Language	25

1. Safety Instructions

-This colony counter has been designed for laboratory usage only.

-Please read this manual carefully before installing or operating this equipment. The instrument shall not be modified in any way. Any modification will void the warranty and may result in potential safety hazard. We are not responsible for any injury or damage caused by modifying the instrument without authorization.

-To avoid the instrument from being burnt or disordered on operation, please use the power adaptor and USB connecting cable supplied by the manufacturer.

-Don't scratch the surface of the Petri dish holder to avoid affecting the illumination of the backlight.

-If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired. The instrument shall not be modified or altered in any way. Any modification or alteration will void the warranty, void the regulatory certifications and create a potential safety hazard.

-We are not responsible for any injury or damage caused by using the instrument for any non-intended purpose, or modifying the instrument by any person who is not authorized.

-Operating environment: Indoor use, 5 °C to 40 °C max. 80% RH max. Main supply voltage fluctuates up to $\pm 10\%$ of the nominal voltage. Overvoltage category I. Pollution degree II.

2. Inspection

2.1 Receiving Inspection

Unpack the equipment carefully and check for any damages which may have arisen during transport. Please contact your supplier for technical support.

Note:

If there is any apparent damage to the system, please do not plug it into the power line.


2.2 Listing of Items

The package includes the following items:

Items in CC100 Colony counter	Qty
Colony counter	1
Magnifier	1
Magnifier holder	1
AC/DC switch adaptor	1
Marking pen (black)	1
Screwdriver	1
User Manual	1
Items in CC100 Software (Optional)	Qty
PC Software USB flash drive	1
USB connecting cable	1
Instruction manual	1

Table 1

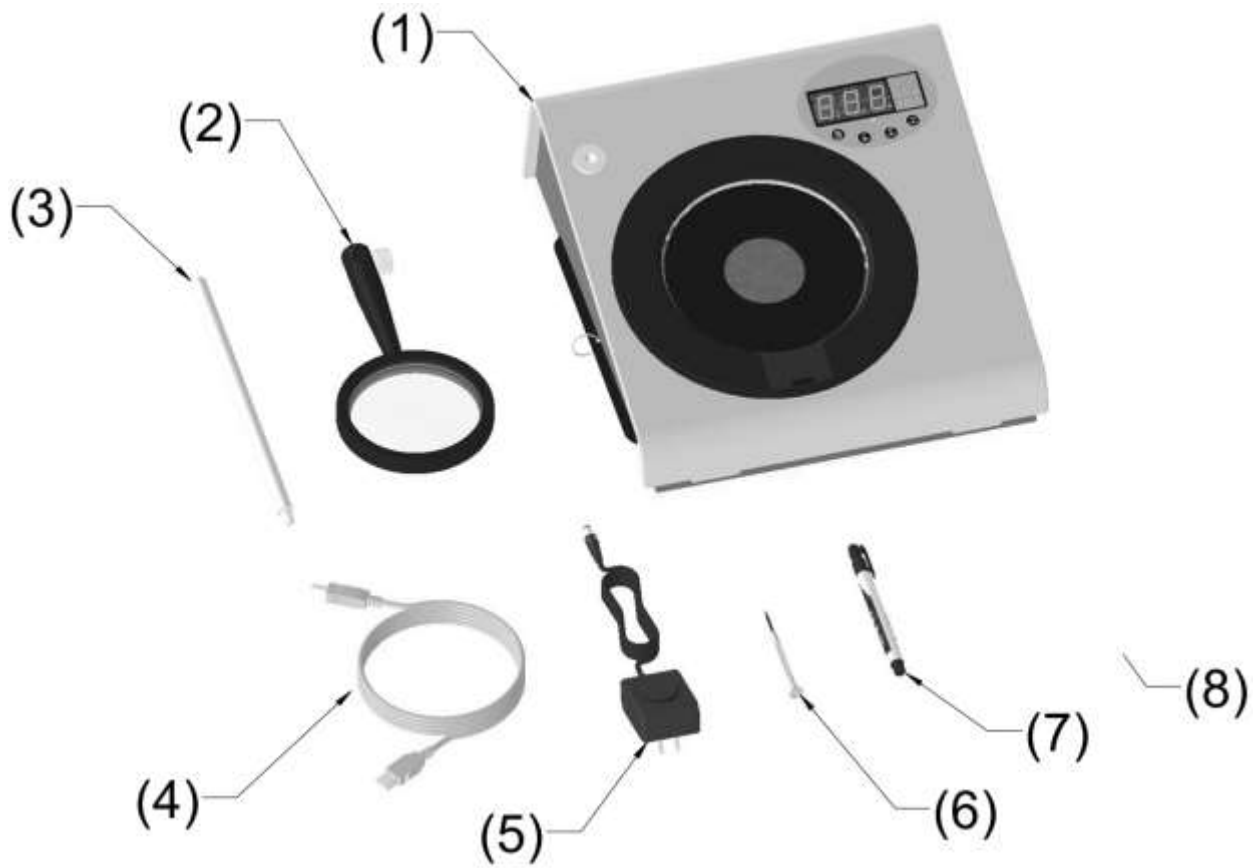
3. Assembly

Item description	Number
Colony counter	1
Magnifier	2
Magnifier holder	3
USB connecting cable (optional)	4
AC/DC switching adaptor	5
Screwdriver	6
Marking pen (black)	7
PC Software USB flash drive (optional)	8

Table 2

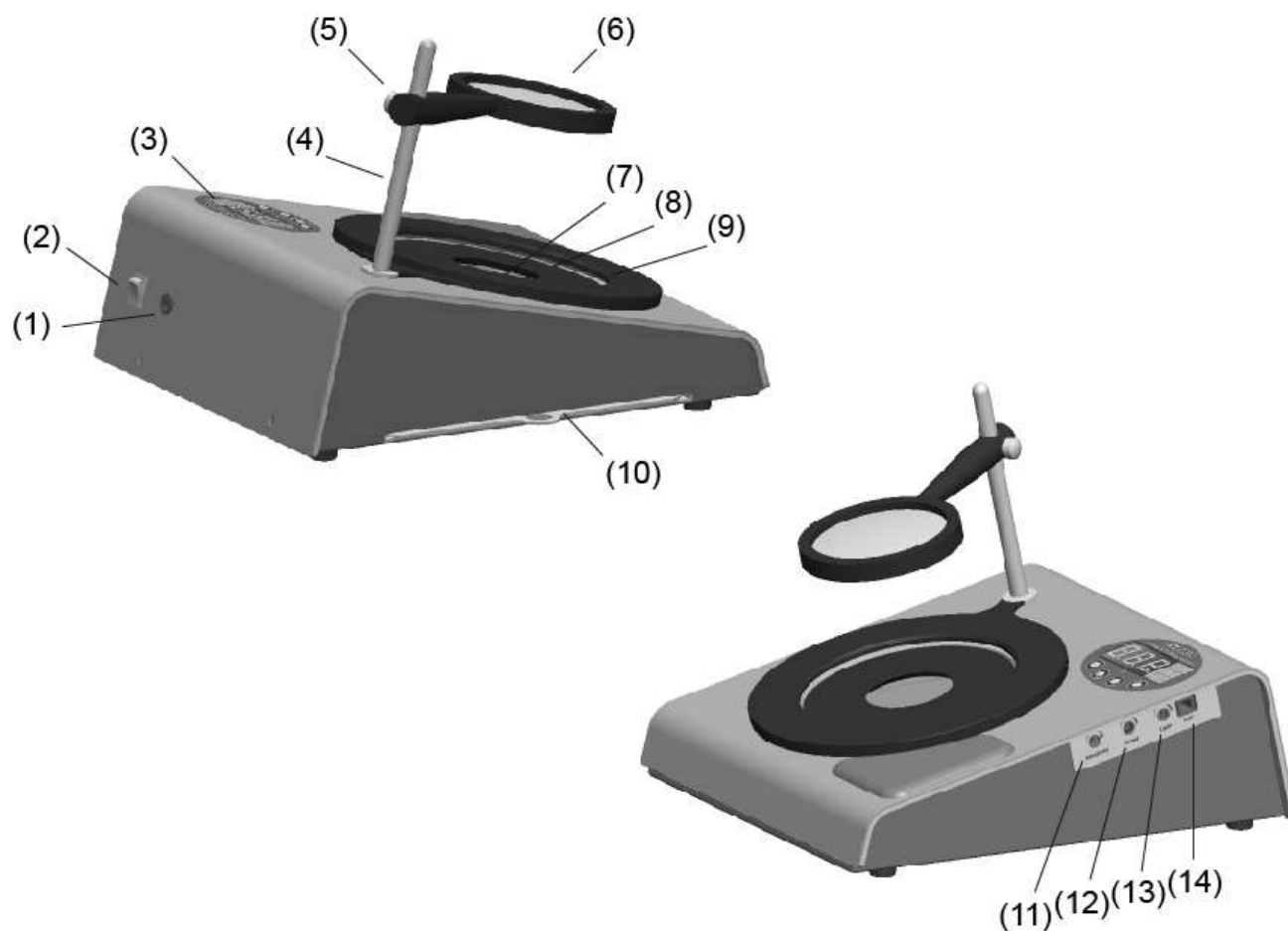
4. Main part diagram

Item description	Number
Power socket	1
Power switch	2
Control panel	3
Magnifier holder	4
Adjustable screw	5
Magnifier	6
Petri dish position	7
Petri dish holder (60 mm)	8
Adjustable petri dish holder (90~150 mm)	9
Background plate	10
Pressure sensitivity regulator	11
Acoustic counting signal regulator	12
Light intensity regulator	13
USB connecting port	14

Table 3

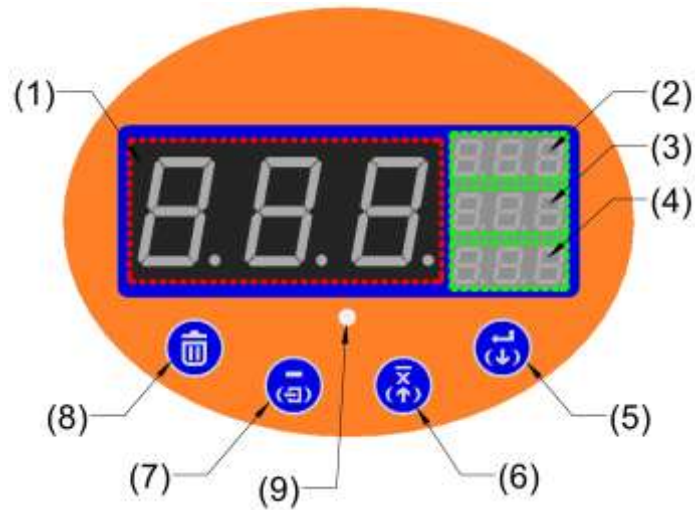
5. Control panel

Item description	Number
Counted value display (LED-A)	1
Storage value display (LED-B)	2
Storage value display (LED-C)	3
Storage value display (LED-D)	4
Save / Next	5
Average / Previous	6
Back / Exit	7
Delete	8
USB connecting indicator (green: connected, red: off)	9

Table 4

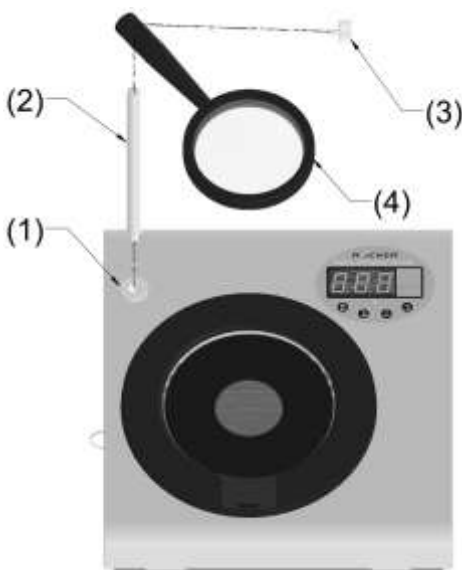6. Installation

Item description	Number
Magnifier position	1
Magnifier holder	2
Adjustable screw	3
Magnifier	4

Table 5

1. Plug the power cord into the main unit and make sure the supplied voltage is within affordable range of power adaptor.
2. Install the magnifier on the main unit by following above diagram and adjust the height of magnifier properly.
3. If you want to transfer the counting data to a computer, please use USB cable provided to connect PC and colony counter and start the operation of software. For the first time use, please install the software by referring to the file of "PC Software Installation and Operation" on USB flash drive.
4. Please use the screwdriver provided to adjust the pressure sensitivity, acoustic volume, illumination required.
5. Change background plate (White/Black) if necessary
6. Please keep the Petri dish position and the magnifier clean to assure good light transmission.

7. Operation

Key and Signal

(1) Normal Mode

	SAVE	Store the count in the memory
	AVERAGE	Average saved counts < display value on (LED-A) >
	BACK	Decrease the value on (LED-A) (alarm with a beep)
	DELETE	Delete the count < alarm with long beep >

(2) View Mode

	EXIT	1). Press and hold this key for 1 second to enter < View Mode > when the display on (LED-A) is "000". 2). Press this key again to exit the <View Mode>
	NEXT	Check the next record
	PREVIOUS	Check the previous record
	DELETE	Delete selected record on display (LED-B/C/D)

(3) Indicator Light

- **Green light:** when the green light is ON means the connection is normal and the data transmission is successful. When light is flashing means the data is under transmission.
- **Red light:** when the red light is ON means the connection is abnormal and data transmission is failing.
- **No light:** it means the PC connection is not detectable.

Operation under Normal Mode

	<p>Place the Petri dish (1) on the dish holder and slide the adjuster (2) properly. Attention: For the Petri dish with 60mm in diameter, you can use “Petri dish holder (60mm)” to fix the Petri dish.</p>
	<p>Ensure the display on LED-A is "000" and no count display on LED-B, LED-C and LED-D before counting. If the count shows up, please press “DELETE” key to clear the memory.</p>
	<p>Mark each colony with the marker pen. Every time a colony is marked, the apparatus will register the count on LED-A with a beep and counter advance.</p>
	<p>Press the “BACK” key if over counting occurs (alarm with a beep)</p>
	<p>After counting, press the “SAVE” key, the saved count will be displayed on LED-B.</p>
	<p>Place the next Petri dish and press <SAVE > key to resume the count. Repeat until all dishes have been counted. The saved counts will display on LED-B, LED-C, LED-D in order.</p>
	<p>After you finish counting, press and hold “EXIT” key for 1 second to enter < View Mode > to check the saved counts. Press “EXIT” key again to be back to <Normal Mode></p>
	<p>Press the <AVERAGE> key to display the average count that calculated from all previously saved counts.</p>

Notice:

- If the counted Petri dishes are over 3 batches, the display will only show the record of latest 3 batches. For the record of previous counts, it will be saved, but not shown on display.
- The main unit can store up to 100 records in memory. It will fail to work if the record is over 100. Please delete the record or activate "AVERAGE" facility, so the main unit can resume working.
- If the main unit connects to PC, the saved counts and average counts will be sent to the software on PC when you press "AVERAGE" key.
- Please refer to the file of "Installation and Operation of Software" on USB flash drive for operation of software.

Operation under View Mode

	<p>Press and hold "EXIT" key for 1 second to enter < View Mode >. Notice: The display LED-A will show the total record numbers before you release the "EXIT" key.</p>
	<p>After releasing the "EXIT" key, the saved counts will display on LED-B/C/D. The selected count will flash and LED-A will show the record number of selected count.</p>
	<p>Press the "NEXT" key to check the next record. It will alarm with a beep when it goes to the last record.</p>
	<p>Press the "PREVIOUS" key to check the previous record. It will alarm with a beep when it goes to the first record.</p>
	<p>Press "DELETE" key to delete flashed record. The alarm will sound with a long beep.</p>
	<p>Press "EXIT" key to be back to <Normal Mode></p>

8. Maintenance

- After every use, please wipe the Petri dish position, the magnifier and surface of colony counter by 75% alcohol for sterilization and to gain good illumination.
- Please keep the instrument in well packed box to avoid dust if is not going to be used plan not to use it for a long time.

9. Troubleshooting

Symptoms	Possible causes and Solution
Device fails to start	1. Check the power supply conform to power specified (DC12V) on main unit 2. Check if the power switch is in the ON position 3. Malfunction of internal parts or internal tubing are loose → contact your distributor for repair
Poor LED lighting	1. Adjust the light intensity with a screwdriver 2. The backlight plate or mainboard is broken → contact your distributor for repair.
Acoustic counting signal is too low	1. Adjust the volume with a screwdriver 2. Mainboard is broken → contact your distributor for repair.
Poor pressure sensitivity	1. Adjust the sensitivity with a screwdriver 2. Mainboard is broken → contact your distributor for repair.
The main unit fails to connect with PC	1. Check if the USB cable is connected to the main unit and computer. 2. Check if the USB driver programming is installed 3. Make sure if the PC software is active. 4. Mainboard is broken → contact your distributor for repair.

10. Software Instructions (Optional)

- This software can only be installed on PC or Laptop running Windows operating system. Please use the USB 2.0 cable supplied with the software package to ensure accurate data transfer.
- Recommended System Requirements:


Operating System	Windows XP SP3 32-bit / XP SP2 64-bit Windows Vista SP2 32-bit / Vista SP2 64-bit Windows 7 SP1 32-bit / 7 SP1 64-bit Windows 8 32-bit / 8 64-bit ; Windows 8.1 32-bit / 8.1 64-bit Windows 10 32-bit / 10 64-bit
CPU	Pentium 4 1.5GHz above
RAM	512MB above (1GB preferred)
Hard Disk Space	Installation 100MB above
Interface	USB Port
Resolution	1024 × 768 Pixels

Connection Status:

- When connected to the computer, the main screen displays  and the GREEN indicator on the main unit lights up.
- When disconnected of the computer, the main screen displays  and the RED indicator on the main unit lights up.

11. Software Installation Guide

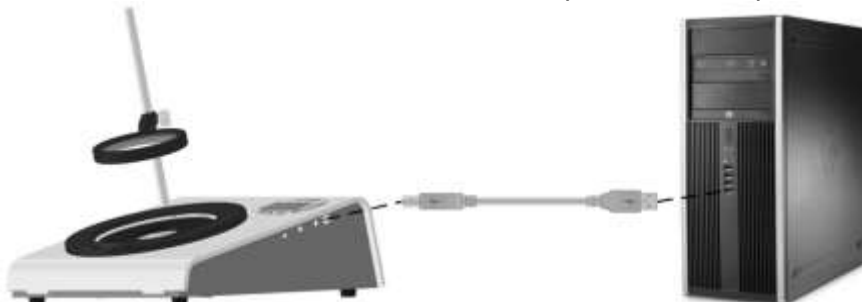- Insert the USB flash disk into one of the USB ports on the computer.

- Execute flash drive : \Galaxy-Setup\Galaxy330Install.exe
- Please click “Allow” or “Yes” if a prompt appears on screen asking for your permission.

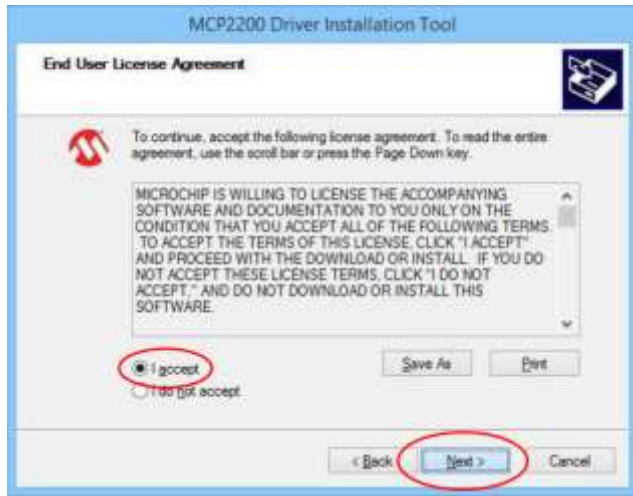
- Click “Setup” to install driver before software installation.
(System auto detect 32- bit or 64 -bit version)


- Note: You must be logged in as an administrator to install this program.

Driver Installation

1. Click “Next” to start installation.
2. Click “I accept” and then click “Next” to begin installation.

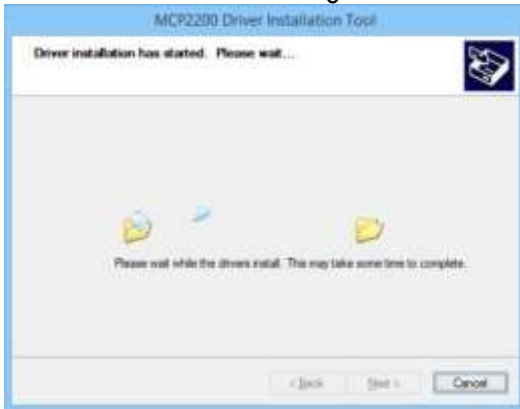3. The driver installation begins and click “Install” to continue installation process.

4. Click "Finish" to complete the installation then begin installing software.

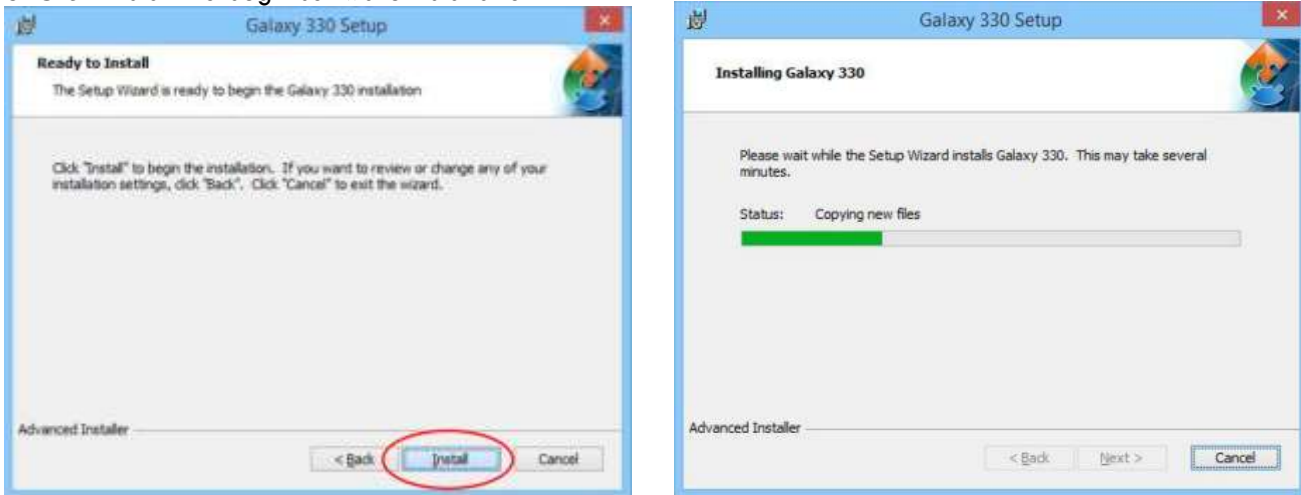
Software Installation

1. Click "Next" to begin software installation.
2. Select installation folder then click "Next" to continue installation process.

3. Click "Install" to begin software installation.

4. Click "Finish" to complete the installation

12. Software operation guide – Data access

(1) Start Program

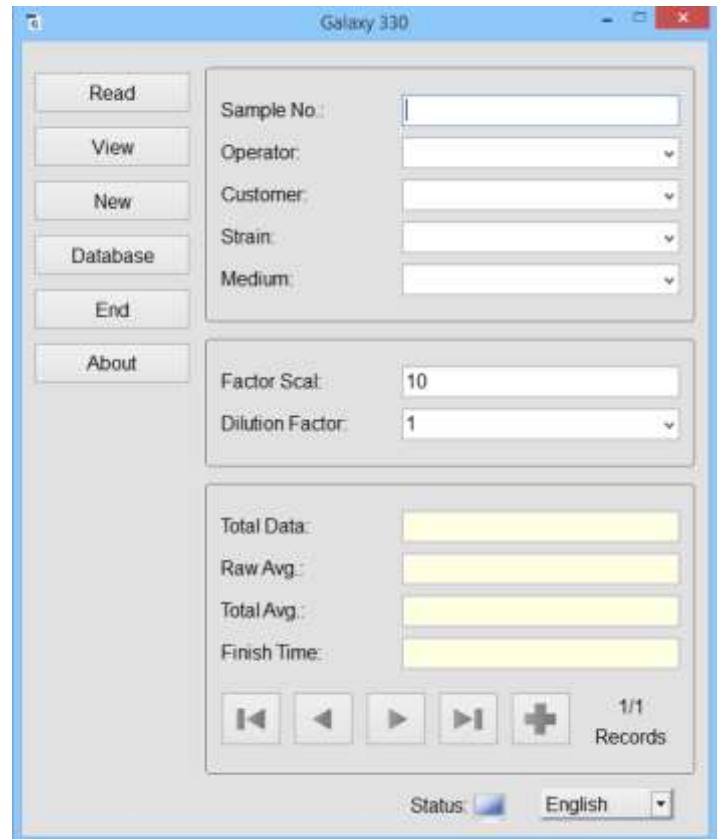
1. Click the Galaxy_330 file icon (as shown below) under the Program Files folder or on the desktop.

2. Execute Galaxy 330 program then the main program screen will be displayed as the right figure.

Note: The main window display may vary depending on the operating system (OS), the user can set the main window properties by performing the following steps:

1. For Windows®, change window properties to Windows classic style.
2. For Windows® XP, click Control Panel → Display Properties → Appearance tab then change the color scheme to Windows classic.
3. For Windows® Vista / 7 / 8 / 8.1 / 10, click Control Panel → Personalization → Theme then change to the Windows Classic theme.


(2) Input Sample data

Sample No: Input Sample Number
Operator: Input Operator Name
Customer: Input Customer Name
Strain: Input Strain Name
Media: Input Media Name
Factor Scal.: Input Dilution Ratio
Dilution Factor: Input Dilution Factor

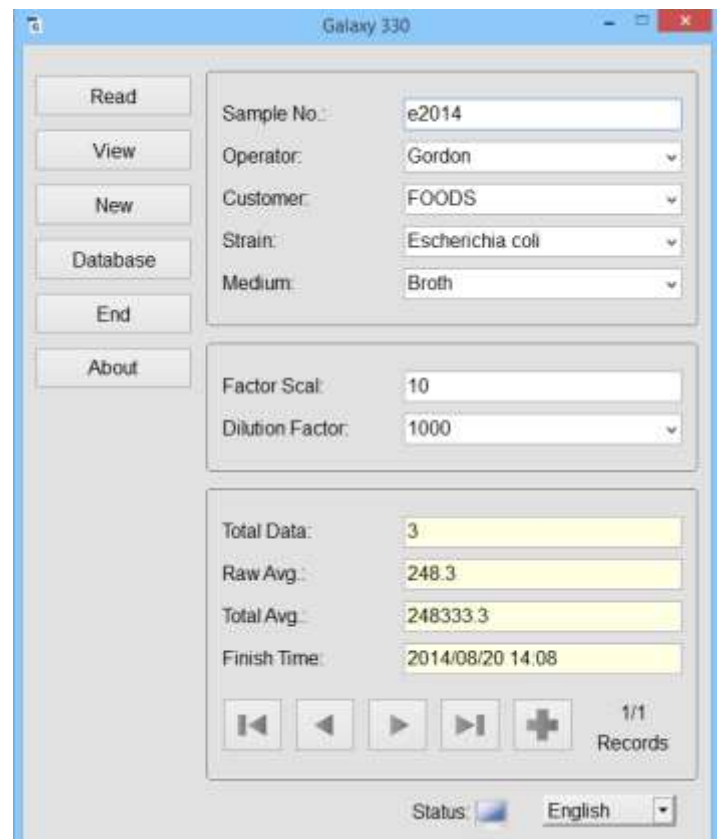(3) Connect to Computer

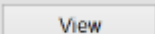
1. Connect to computer with USB cable


- Switch on Colony Counter then the Status indicator will display connection signal.
- When connected, the status indicator will display  and Colony Counter indicator will display "Green".
- When disconnected, the status indicator will display  and Colony Counter indicator will display "Red".

Note: When the USB is connected, Windows® will automatically detect the hardware and perform driver, please be patient while in process.

(4) Reading and Querying Data

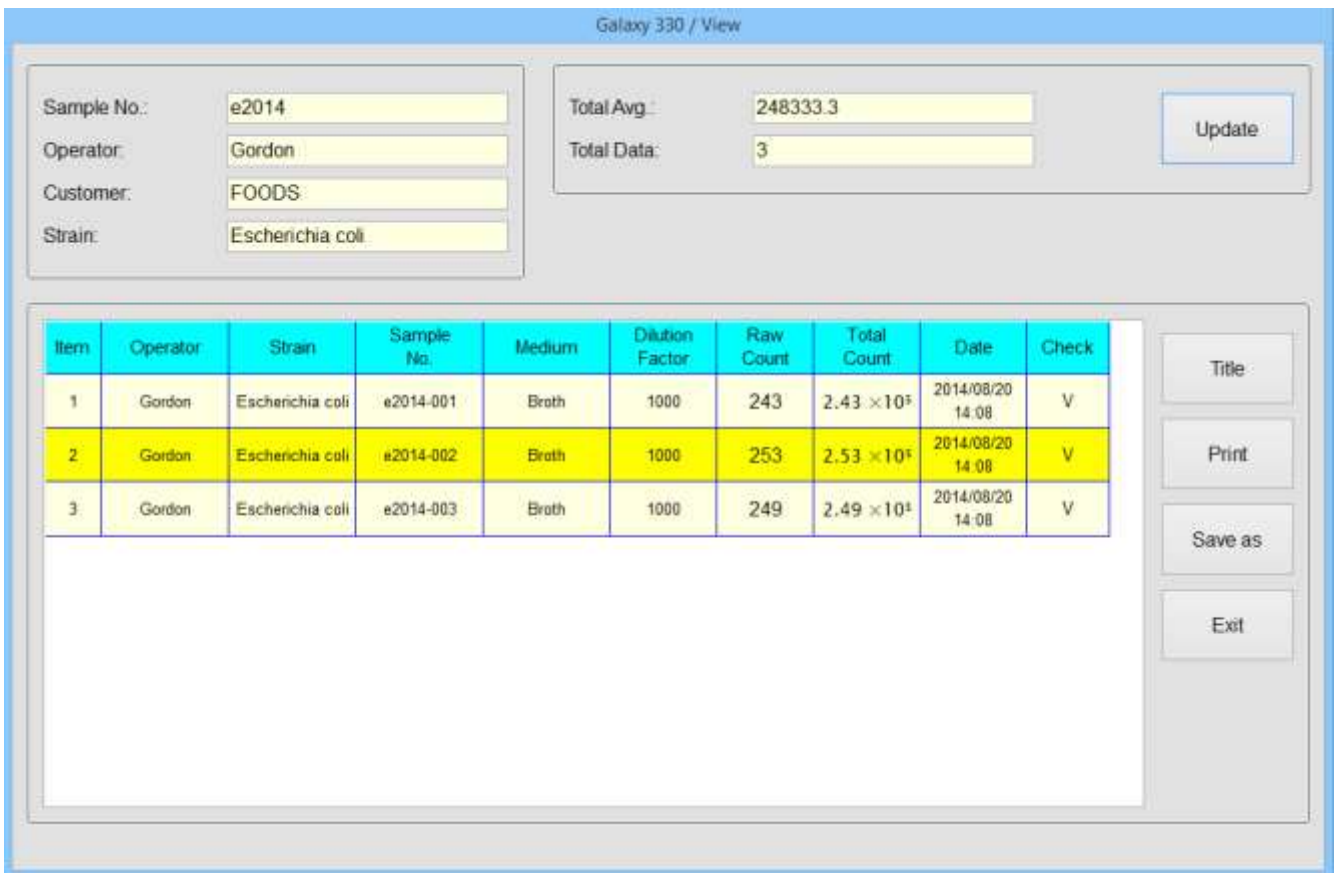- After counting bacterial colonies in the same sample and same dilution factor on each plate at Colony Counter, press "Average Button"  then each average count value will be automatically transmitted to the computer.
- Press  the main program screen and query all data received.

Item	Operator	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date	Check
1	Gordon	Escherichia coli	e2014-001	Broth	1000	243	2.43×10^5	2014/08/20 14:08	V
2	Gordon	Escherichia coli	e2014-002	Broth	1000	253	2.53×10^5	2014/08/20 14:08	V
3	Gordon	Escherichia coli	e2014-003	Broth	1000	249	2.49×10^5	2014/08/20 14:08	V

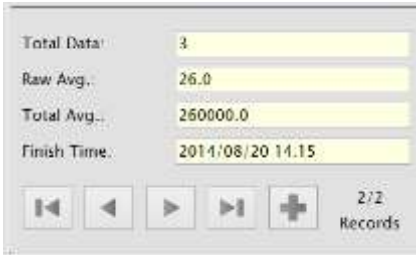
- Press "EXIT" to return to the main program screen.

Note: a set of data can contain max. 100 records and the number from 001 to 999 will be automatically added on counting data for different plate.


Note: in case the field cannot display all complete information, "... " will appear at the end of the field, then move mouse to the field, all complete information will be displayed after one second.

(5) Reading a new data

1. After completion of counting the following plates of bacterial colonies in the same sample and dilution factor at Colony Counter, press "Average Button"  then the next average counting data will be automatically transmitted to the computer again.
2. When turning into 2/2 Records, press each of the four keys below to switch display on counting data in different dilution factors up to 10 sets of data.

	To the first data	
	To the previous data	
	To the next data	
	To the last data	

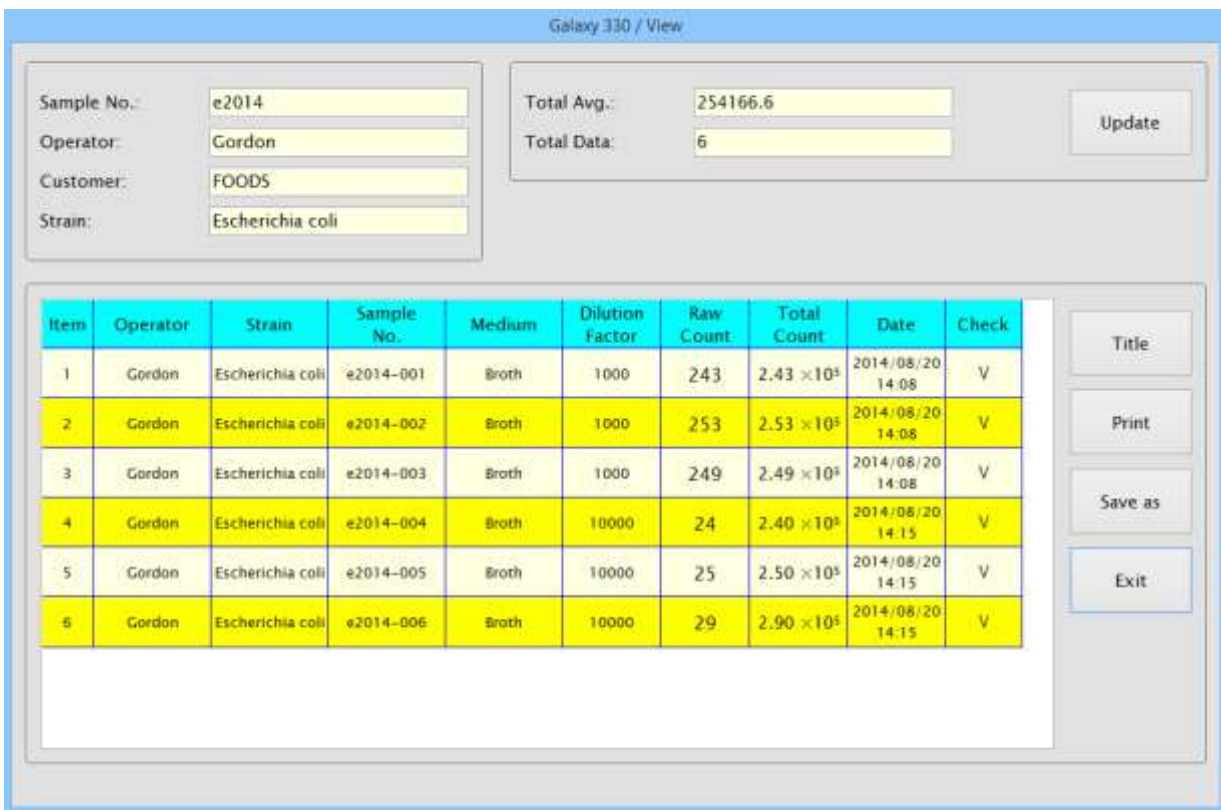3. Press  the main program screen to query all data received.

Galaxy 330 / View

Sample No.: e2014
 Operator: Gordon
 Customer: FOODS
 Strain: Escherichia coli

Total Avg.: 254166.6
 Total Data: 6

Update

Item	Operator	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date	Check
1	Gordon	Escherichia coli	e2014-001	Broth	1000	243	2.43 × 10 ⁵	2014/08/20 14:08	V
2	Gordon	Escherichia coli	e2014-002	Broth	1000	253	2.53 × 10 ⁵	2014/08/20 14:08	V
3	Gordon	Escherichia coli	e2014-003	Broth	1000	249	2.49 × 10 ⁵	2014/08/20 14:08	V
4	Gordon	Escherichia coli	e2014-004	Broth	10000	24	2.40 × 10 ⁵	2014/08/20 14:15	V
5	Gordon	Escherichia coli	e2014-005	Broth	10000	25	2.50 × 10 ⁵	2014/08/20 14:15	V
6	Gordon	Escherichia coli	e2014-006	Broth	10000	29	2.90 × 10 ⁵	2014/08/20 14:15	V

Title
 Print
 Save as
 Exit

Press "EXIT" to return to the main program screen.

(6) Viewing Data

Please repeat the previous steps for counting the same sample in different dilution factors. After completion of counting all different dilution factors for the same sample, press "View" to view all counting data, delete data, print data or save data as Excel file. Press "Exit" to return to the main program screen.

<<Please refer to respective instruction on how to delete, print or save data below>>

Note: in case the field cannot display all the information, "... " will appear at the end of the field, then move mouse to the field, all complete information will be displayed after one second.

(7) Completing the count

After completing the counts of the previous sample, please press  and follow the above (1) ~ (6) steps to count dilution factors for the next sample.

Note:

- a. After pressing the "New" key, a confirmation screen will appear. Press confirmation key to store all received data in database and then press to view the set of data just saved;
- b. This software can store up to max. 100,000 sets of data, once the number of the stored data exceeds the max. number, it will automatically delete the earliest data and then store the newly created data.

13. Software operation guide – Data deletion

1- Enter View Screen

During connection process, press to query all received data in total 10 fields.

Note: in case the field cannot display all complete information, "..." will appear at the end of the field, then move mouse to the field, all complete information will be displayed after one second.

Item	Operator	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date	Check
1	Gordon	Escherichia coli	e2014-001	Synthetic me...	1000	243	2.43 ×10 ⁵	2014/08/20 14:08	V
2	Gordon	Escherichia coli	e2014-002	Synthetic me...	1000	253	2.53 ×10 ⁵	2014/08/20 14:08	V
3	Gordon	Escherichia coli	e2014-003	Synthetic me...	1000	249	2.49 ×10 ⁵	2014/08/20 14:08	V
4	Gordon	Escherichia coli	e2014-004	Synthetic me...	10000	24	2.40 ×10 ⁵	2014/08/20 14:15	V
5	Gordon	Escherichia coli	e2014-005	Synthetic me...	10000	25	2.50 ×10 ⁵	2014/08/20 14:15	V
6	Gordon	Escherichia coli	e2014-006	Synthetic me...	10000	29	2.90 ×10 ⁵	2014/08/20 14:15	V

Item	Data Number
Operator	Operator Name
Strain	Strain Name
Sample No	Counted Sample Number
Media	Media Name
Dilution Factor	Dilution Factor for a concentrated sample
Raw Count	Colony count number
Total Count	The number of colonies after conversion
Date	The data-receiving date
Check	Check if the data is valid

2- Select the data you want to delete

If any incorrect count data is found, move mouse cursor to the Check field.

Item	Operator	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date	Check
1	Gordon	Escherichia coli	e2014-001	Synthetic me...	1000	243	2.43 ×10 ⁵	2014/08/20 14:08	V
2	Gordon	Escherichia coli	e2014-002	Synthetic me...	1000	253	2.53 ×10 ⁵	2014/08/20 14:08	V
3	Gordon	Escherichia coli	e2014-003	Synthetic me...	1000	249	2.49 ×10 ⁵	2014/08/20 14:08	V
4	Gordon	Escherichia coli	e2014-004	Synthetic me...	10000	24	2.40 ×10 ⁵	2014/08/20 14:15	V
5	Gordon	Escherichia coli	e2014-005	Synthetic me...	10000	25	2.50 ×10 ⁵	2014/08/20 14:15	V
6	Gordon	Escherichia coli	e2014-006	Synthetic me...	10000	29	2.90 ×10 ⁵	2014/08/20 14:15	V

Click the left mouse button to "V" sign to make the field become blank

Item	Operator	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date	Check
1	Gordon	Escherichia coli	e2014-001	Synthetic me...	1000	243	2.43 × 10 ⁵	2014/08/20 14:08	V
2	Gordon	Escherichia coli	e2014-002	Synthetic me...	1000	253	2.53 × 10 ⁵	2014/08/20 14:08	
3	Gordon	Escherichia coli	e2014-003	Synthetic me...	1000	249	2.49 × 10 ⁵	2014/08/20 14:08	V
4	Gordon	Escherichia coli	e2014-004	Synthetic me...	10000	24	2.40 × 10 ⁵	2014/08/20 14:15	V
5	Gordon	Escherichia coli	e2014-005	Synthetic me...	10000	25	2.50 × 10 ⁵	2014/08/20 14:15	V
6	Gordon	Escherichia coli	e2014-006	Synthetic me...	10000	29	2.90 × 10 ⁵	2014/08/20 14:15	

Note: The blank field indicates the data will be deleted and re-click on the blank field will show the following screen. Click on the unwanted data and press "Update" to confirm the deletion.

Sample No.:	<input type="text" value="e2014"/>	Total Avg.:	<input type="text" value="245500.0"/>	<div style="border: 2px solid red; padding: 5px; display: inline-block;">Update</div>
Operator:	<input type="text" value="Gordon"/>	Total Data:	<input type="text" value="4"/>	
Customer:	<input type="text" value="FOODS"/>			
Strain:	<input type="text" value="Escherichia coli"/>			

Press "Yes" to reconfirm the data deletion. (unable to be recovered after deletion)

Galaxy 330 / Warning

Change Data!! Yes/No

Yes

No

Shown as below after data deletion.

Galaxy 330 / View

Sample No.:	<input type="text" value="e2014"/>	Total Avg.:	<input type="text" value="245500.0"/>	<div style="border: 1px solid gray; padding: 5px; display: inline-block;">Update</div>
Operator:	<input type="text" value="Gordon"/>	Total Data:	<input type="text" value="4"/>	
Customer:	<input type="text" value="FOODS"/>			
Strain:	<input type="text" value="Escherichia coli"/>			

Item	Operator	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date	Check
1	Gordon	Escherichia coli	e2014-001	Synthetic me...	1000	243	2.43 × 10 ⁵	2014/08/20 14:08	V
2	Gordon	Escherichia coli	e2014-003	Synthetic me...	1000	249	2.49 × 10 ⁵	2014/08/20 14:08	V
3	Gordon	Escherichia coli	e2014-004	Synthetic me...	10000	24	2.40 × 10 ⁵	2014/08/20 14:15	V
4	Gordon	Escherichia coli	e2014-005	Synthetic me...	10000	25	2.50 × 10 ⁵	2014/08/20 14:15	V

Title

Print

Save as

Exit

(3) After completion of data deletion and press "Exit" to return to the main program screen.

14. Software operation guide – Data Query

-Enter View Screen

Press "Database" in the main program screen to enter view screen.

Galaxy 330 / Database

Sample No.:

Operator:

Customer:

Strain:

Start Date: 2014/08/20

End Date: 2014/08/20

Total Data:

Total Avg.:

Search

Clear

Item	Operator	Customer	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date
------	----------	----------	--------	------------	--------	-----------------	-----------	-------------	------

Title

Print

Save as

Exit

The six fields below can be set as query conditions.

Sample No.
Operator
Customer
Strain
Start Date
End Date

Note: The query condition can accept English capital, small letters or incomplete characters.

-Start Search

Press "Search" after inputting the query condition then all queried data will appear.

Press "Clear" to clear all displayed data before re-query.

Galaxy 330 / Database

Sample No.:

Operator: Gordon

Customer:

Strain:

Start Date: 2014/08/20

End Date: 2014/08/20

Total Data: 4

Total Avg.: 245500.0

Search

Clear

Item	Operator	Customer	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date
1	Gordon	FOODS	Escherichia...	e2014-001	Synthetic me...	1000	243	2.43×10^5	2014/08/20 14:08
2	Gordon	FOODS	Escherichia...	e2014-003	Synthetic me...	1000	249	2.49×10^5	2014/08/20 14:08
3	Gordon	FOODS	Escherichia...	e2014-004	Synthetic me...	10000	24	2.40×10^5	2014/08/20 14:15
4	Gordon	FOODS	Escherichia...	e2014-005	Synthetic me...	10000	25	2.50×10^5	2014/08/20 14:15

Title

Print

Save as

Exit

Note: This software can display up to 1000 sets of data. Please narrow your search to avoid excessive data.

-Search Result

The search result will be displayed as above in total 10 fields incl. Item No., Operator, Customer, Strain, Sample No, Media, Dilution Factor, Raw Count, Total Count and Date.

Note: in case the field cannot display all complete information, "..." will appear at the end of the field, then move mouse to the field, all complete information will be displayed after one second.

15. Software operation guide – Data Print

(1) Search the data for printing

Search the data for printing by pressing “View” or “Database” in the main program screen and a sub-screen will appear. Press “Print” and set the printer in portrait, A4 format then the data list will be printed out.

The screenshot shows the 'Galaxy 330 / View' software interface. It features a search form at the top with fields for Sample No., Operator, Customer, and Strain. Below the search form is a table with 10 columns: Item, Operator, Strain, Sample No., Medium, Dilution Factor, Raw Count, Total Count, Date, and Check. The table contains 4 rows of data. To the right of the table is a vertical menu with buttons for Title, Print, Save as, and Exit. The 'Print' button is highlighted with a red box.

Item	Operator	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date	Check
1	Gordon	Escherichia coli	e2014-001	Synthetic me...	1000	243	2.43 × 10 ⁵	2014/08/20 14:08	V
2	Gordon	Escherichia coli	e2014-003	Synthetic me...	1000	249	2.49 × 10 ⁵	2014/08/20 14:08	V
3	Gordon	Escherichia coli	e2014-004	Synthetic me...	10000	24	2.40 × 10 ⁵	2014/08/20 14:15	V
4	Gordon	Escherichia coli	e2014-005	Synthetic me...	10000	25	2.50 × 10 ⁵	2014/08/20 14:15	V

The screenshot shows the 'Print' dialog box. It has a 'Printer' section with a dropdown menu set to 'Microsoft XPS Document Writer' and a 'Properties...' button. Below this, it shows the printer's status as 'Ready', type as 'Microsoft XPS Document Writer v4', and location as 'PORTPROMPT:'. There is a 'Print range' section with radio buttons for 'All', 'Pages' (with 'from:' and 'to:' fields), and 'Selection'. The 'Copies' section has a 'Number of copies' spinner set to '1' and a 'Collate' checkbox. At the bottom, there are 'OK' and 'Cancel' buttons.

COLONY COUNTING REPORT

Item	Operator	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date
1	Gordon	Escherichia coli	e2014-001	Synthetic me...	1000	243	2.43 × 10 ⁵	2014/08/20 14:08
2	Gordon	Escherichia coli	e2014-003	Synthetic me...	1000	249	2.49 × 10 ⁵	2014/08/20 14:08
3	Gordon	Escherichia coli	e2014-004	Synthetic me...	10000	24	2.40 × 10 ⁵	2014/08/20 14:15
4	Gordon	Escherichia coli	e2014-005	Synthetic me...	10000	25	2.50 × 10 ⁵	2014/08/20 14:15
Total data: 4					Average: 2.455 × 10 ⁵			

Print date: 2014/08/20 15:46

1

Note: The printing format is designed for the portrait, A4 paper only so please don't use another size. To print the other paper formats. Please save the data as Excel file and edit the file with Excel before printing.

(2) Printing the report with title

Press "Title" in the sub-screen from pressing "View" or "Database" to print the report with title.

Fill in all required information after the following form appears.

Galaxy 330 / Title

Company:

Department: Sales Department

Tel:

Fax:

Address:

E-mail:

Operator: Gordon

Approval: Tim

Print Title

Submit

Cancel

Load

Save

Make sure that "Print Title" has been ticked after filling in all the title information.

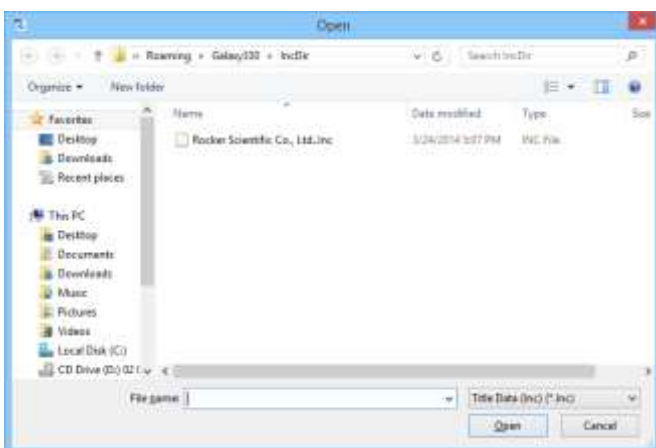
Note: if "Print Title" is not ticked then the title won't be printed out.

(4) Saving Data

Press "Save" to save all title information.

All saved title information can then be re-loaded back into the program by pressing "Load".

(5) Print the report with title

Ensure all title information are correct and press "Submit" to return to the previous screen. After pressing "Print" and selecting the printer, the report with title can be printed out.

Note:

- Press "Cancel" to abort any amended title information and all previous title information will keep unchanged.
- Uncheck "Print Title" if the title information is not required for the report.

COLONY COUNTING REPORT

Company:

Department: Sales Department

Address:

Tel:

Fax:

E-mail:

Operator: Gordon

Approval: Tim


Item	Operator	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date
1	Gordon	Escherichia coli	e2014-001	Synthetic me...	1000	243	2.43 × 10 ⁶	2014/08/20 14:08
2	Gordon	Escherichia coli	e2014-003	Synthetic me...	1000	249	2.49 × 10 ⁶	2014/08/20 14:08
3	Gordon	Escherichia coli	e2014-004	Synthetic me...	10000	24	2.40 × 10 ⁶	2014/08/20 14:13
4	Gordon	Escherichia coli	e2014-005	Synthetic me...	10000	25	2.50 × 10 ⁶	2014/08/20 14:13
Total data: 4					Average: 2.455 × 10 ⁶			

Print date: 2014/08/20 15:48

16. Software operation guide – Data Saving

(1) Search the data to be saved

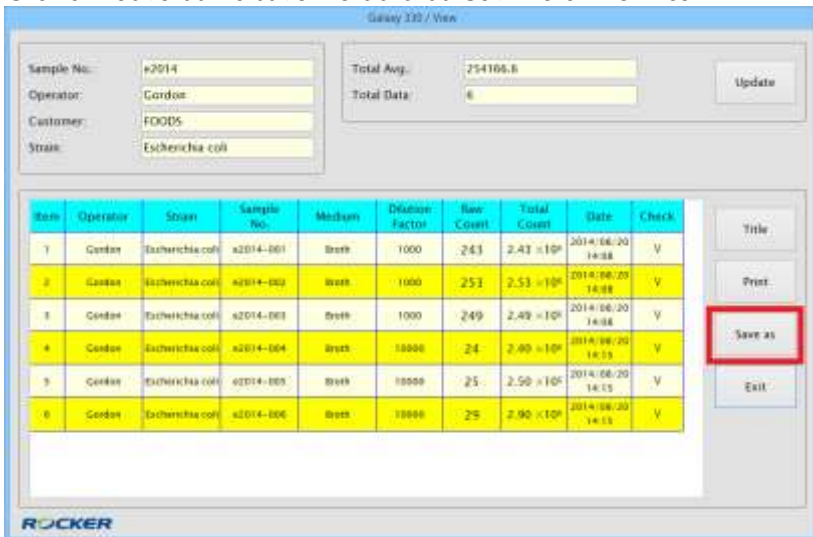Search the data to be saved through the sub-screen from pressing "View" or "Database".
Click on "Save as" to save the data as CSV file of the Excel.

(2) Open .CSV File with Excel

After opening Excel, select open the old file (.csv file).

Item	Operator	Customer	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date
1	Gordon	FOGDS	Escherichia coli	e2014-001	Synthetic media	1000	243	2.43E+05	2014/8/20 14:08
2	Gordon	FOGDS	Escherichia coli	e2014-003	Synthetic media	1000	249	2.49E+05	2014/8/20 14:08
3	Gordon	FOGDS	Escherichia coli	e2014-004	Synthetic media	10000	24	2.40E+05	2014/8/20 14:15
4	Gordon	FOGDS	Escherichia coli	e2014-005	Synthetic media	10000	25	2.50E+05	2014/8/20 14:15

After opening the .csv file, all the data can be tidied as desired.

Note:

- a. Whether ticking “Print Title” or not, all existing title information will be included in the .CSV file. However, these title information can be edited or deleted as requested.
- b. The data of total count will be expressed in scientific notation.

17. Software operation guide – Change Language

CHANGE LANGUAGE:

Move mouse cursor to the right button at the main program screen to switch English / Traditional Chinese/Simple Chinese lanaguage interfaces:

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The screenshot shows the 'Galaxy 330 / Database' window. It features a search form with the following fields: Sample No., Operator (dropdown), Customer (dropdown), Strain (dropdown), Start Date (calendar), End Date (calendar), Total Data (text), and Total Avg. (text). There are 'Search' and 'Clear' buttons. Below the search form is a data table with the following columns: Item, Operator, Customer, Strain, Sample No., Medium, Dilution Factor, Raw Count, Total Count, and Date. To the right of the table are buttons for 'Title', 'Print', 'Save as', and 'Exit'.

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The screenshot shows the 'Galaxy 330 / Database' window after a search. The search criteria are: Sample No. (empty), Operator: Gordon, Customer (empty), Strain (empty), Start Date: 2014/08/20, End Date: 2014/08/20, Total Data: 4, and Total Avg.: 245500.0. The 'Search' and 'Clear' buttons are highlighted with a red box. The data table below shows the following results:

Item	Operator	Customer	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date
1	Gordon	FOODS	Escherichia...	e2014-001	Synthetic me...	1000	243	2.43 × 10 ⁵	2014/08/20 14:08
2	Gordon	FOODS	Escherichia...	e2014-003	Synthetic me...	1000	249	2.49 × 10 ⁵	2014/08/20 14:08
3	Gordon	FOODS	Escherichia...	e2014-004	Synthetic me...	10000	24	2.40 × 10 ⁵	2014/08/20 14:15
4	Gordon	FOODS	Escherichia...	e2014-005	Synthetic me...	10000	25	2.50 × 10 ⁴	2014/08/20 14:15

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The screenshot shows the 'Galaxy 330 / View' software interface. At the top, there are input fields for 'Sample No.' (e2014), 'Operator' (Gordon), 'Customer' (FOODS), and 'Strain' (Escherichia coli). To the right, there are summary fields for 'Total Avg.' (245500.0) and 'Total Data' (4), with an 'Update' button. Below this is a table with the following data:

Item	Operator	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date	Check
1	Gordon	Escherichia coli	e2014-001	Synthetic me...	1000	243	2.43 × 10 ⁵	2014/08/20 14:08	V
2	Gordon	Escherichia coli	e2014-003	Synthetic me...	1000	249	2.49 × 10 ⁵	2014/08/20 14:08	V
3	Gordon	Escherichia coli	e2014-004	Synthetic me...	10000	24	2.40 × 10 ⁵	2014/08/20 14:15	V
4	Gordon	Escherichia coli	e2014-005	Synthetic me...	10000	25	2.50 × 10 ⁵	2014/08/20 14:15	V

On the right side of the table, there is a vertical menu with buttons: 'Title', 'Print' (highlighted with a red box), 'Save as', and 'Exit'. An 'Update' button is also present in the top right corner.

The screenshot shows the 'Print' dialog box. It has a title bar with 'Print' and a close button. The dialog is divided into several sections:

- Printer:** Name: Microsoft XPS Document Writer (with a dropdown arrow and 'Properties...' button). Status: Ready. Type: Microsoft XPS Document Writer v4. Where: PORTPROMPT. Comment: (empty).
- Print range:** Radio buttons for 'All' (selected), 'Pages' (with 'from:' and 'to:' input fields), and 'Selection'.
- Copies:** 'Number of copies:' set to 1 (with a spinner). A 'Collate' checkbox is present and unchecked. Below this are three document icons labeled 1, 2, and 3.
- Buttons:** 'OK' and 'Cancel' buttons at the bottom.

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The screenshot shows the Galaxy 530 software interface. At the top, there are input fields for Sample No. (e2014), Operator (Gordon), Customer (FOODS), and Strain (Escherichia coli). To the right, there are summary statistics: Total Avg.: 245500.0 and Total Data: 4, with an Update button. Below this is a data table with columns: Item, Operator, Strain, Sample No., Medium, Dilution Factor, Raw Count, Total Count, Date, and Check. The table contains 4 rows of data. To the right of the table is a sidebar with buttons: Title (highlighted with a red box), Print, Save as, and Exit. The Rocker logo is visible in the bottom left corner.

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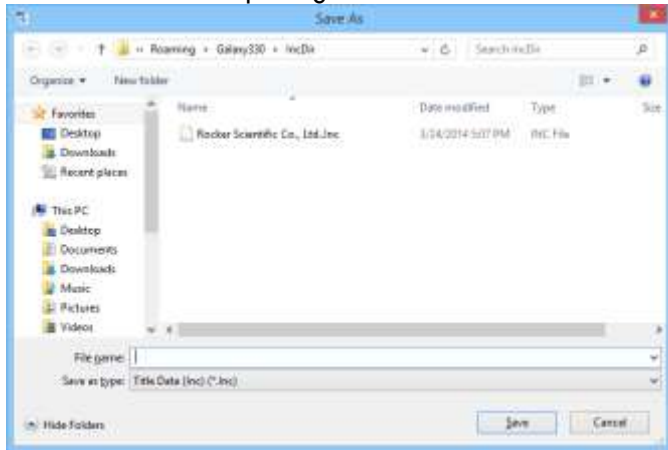
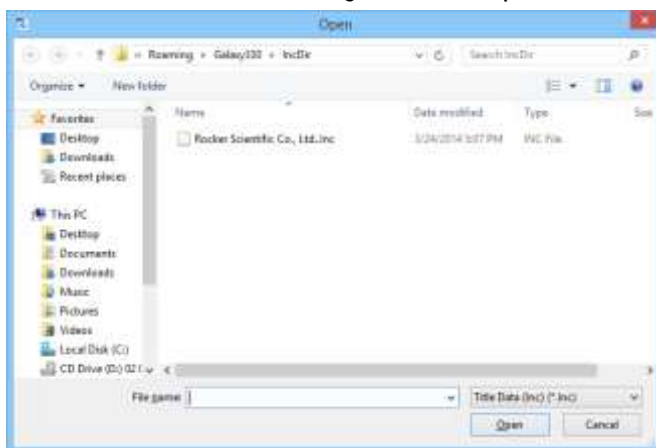
The screenshot shows the 'Colony Counting Software' interface. On the left, there are input fields for Sample No. (e2014), Operator (Gordon), Customer (FOODS), and Strain (Escherichia coli). In the center, there are summary statistics: Total Avg. (254106.6) and Total Data (6). On the right, there are buttons for 'Update', 'Print', 'Save as', and 'Exit'. Below these is a table with columns: Item, Operator, Strain, Sample No., Medium, Dilution Factor, Raw Count, Total Count, Date, and Check. The table contains 6 rows of data. To the right of the software window, a 'Save As' dialog box is open, showing the file name '20140820' and the save type 'Excel 2012 file (*.xlsx)'. The 'Save as' button in the software interface is highlighted with a red box.

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The screenshot shows the 'Galaxy 330 / Database' window. It features a search criteria section at the top with fields for Sample No., Operator, Customer, Strain, Start Date, End Date, Total Data, and Total Avg. Below this is a large table with columns: Item, Operator, Customer, Strain, Sample No., Medium, Dilution Factor, Raw Count, Total Count, and Date. To the right of the table are buttons for Title, Print, Save as, and Exit.

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This screenshot shows the same software interface as above, but with search results displayed in the table. The 'Search' and 'Clear' buttons are highlighted with a red box. The table contains four rows of data.

Item	Operator	Customer	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date
1	Gordon	FOODS	Escherichia ...	e2014-001	Synthetic me...	1000	243	2.43 × 10 ⁵	2014/08/20 14:08
2	Gordon	FOODS	Escherichia ...	e2014-003	Synthetic me...	1000	249	2.49 × 10 ⁵	2014/08/20 14:08
3	Gordon	FOODS	Escherichia ...	e2014-004	Synthetic me...	10000	24	2.40 × 10 ⁵	2014/08/20 14:15
4	Gordon	FOODS	Escherichia ...	e2014-005	Synthetic me...	10000	25	2.50 × 10 ⁴	2014/08/20 14:15

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The screenshot shows the 'Galaxy 330 / View' software interface. At the top, there are input fields for 'Sample No.' (e2014), 'Operator' (Gordon), 'Customer' (FOODS), and 'Strain' (Escherichia coli). To the right, there are summary statistics: 'Total Avg.: 245500.0' and 'Total Data: 4', with an 'Update' button. Below this is a table with 10 columns: Item, Operator, Strain, Sample No., Medium, Dilution Factor, Raw Count, Total Count, Date, and Check. The table contains 4 rows of data. To the right of the table is a vertical menu with buttons for 'Title', 'Print' (highlighted with a red box), 'Save as', and 'Exit'.

Item	Operator	Strain	Sample No.	Medium	Dilution Factor	Raw Count	Total Count	Date	Check
1	Gordon	Escherichia coli	e2014-001	Synthetic me...	1000	243	2.43 × 10 ⁵	2014/08/20 14:08	V
2	Gordon	Escherichia coli	e2014-003	Synthetic me...	1000	249	2.49 × 10 ⁵	2014/08/20 14:08	V
3	Gordon	Escherichia coli	e2014-004	Synthetic me...	10000	24	2.40 × 10 ⁵	2014/08/20 14:15	V
4	Gordon	Escherichia coli	e2014-005	Synthetic me...	10000	25	2.50 × 10 ⁵	2014/08/20 14:15	V

The screenshot shows the Windows 'Print' dialog box. The 'Printer' section shows 'Microsoft XPS Document Writer' as the selected printer, with a 'Properties...' button. The status is 'Ready' and the type is 'Microsoft XPS Document Writer v4'. The 'Print range' section has 'All' selected. The 'Copies' section shows 'Number of copies: 1' and a 'Collate' checkbox. There are 'OK' and 'Cancel' buttons at the bottom.

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The screenshot shows the Galaxy 330 software interface. At the top, there are input fields for Sample No. (e2014), Operator (Gordon), Customer (FOODS), and Strain (Escherichia coli). To the right, there are summary fields for Total Avg. (245500.0) and Total Data (4), with an Update button. Below this is a table with columns: Item, Operator, Strain, Sample No., Medium, Dilution Factor, Raw Count, Total Count, Date, and Check. The table contains four rows of data. To the right of the table is a vertical menu with buttons: 'Titre' (highlighted with a red box), 'Print', 'Save as', and 'Exit'. The Rocker logo is visible in the bottom left corner.

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