ENGLISH - Face shield USER GUIDE

EU – Type certification issued by:

2754: ALIENOR CERTIFICATION

(notified body: 2754) -ZA du Sanital

21 Rue Albert Einstein,**86100 Chatellerault – France****USER INSTRUCTIONS AS PER EU 2016/425****WARNING:**

- The model face protector does not protect against molten metal, hot solids, electric hazard or infrared radiation.

- Materials which might be in contact with the user skin, could cause allergic reactions to susceptible individuals.

- If the face protector is worn over standard ophthalmic spectacles, could transmit impacts of highspeed particles, creating a hazard to the user.

REQUIREMENT: Face protector complies with the requirements of the Regulation EU 2016/425 by means of the European Standard EN 166:2001 "Personal eye protection. Specifications".**PERFORMANCE:** Personal eye protection. Specifications in accordance with EN 166:2001 EU type Examination.**APPLICATION:** When the face protector is worn properly and is given an adequate care, it provides an excellent protection by preventing harmful particles such as dust, grass, flying objects that comes from crafting, cutting, motor saws etc.**MARKINGS:** The appropriate marking can be found on or under the headgear and visor of each model.**VISOR:** CH 1 B  **HEADGEAR:** CH 166 B 3  23865 Taipei Taiwan

Abbreviation as:

CH: Manufacturer Code

B: Medium Mechanical

166: EN Standard

Strength3: Splash Resistant

1: Optical Class (Continuous use) CE: CE Mark

23865: Manufacturer Postal Code

Taipei, Taiwan: Manufacturer City, Country

Optical Classes

Optical class 1 is for work with particularly high vision requirements for permanent use

Optical class 2 is for work with average vision requirements

Optical class 3 is for rough work without any special vision requirements (only intended for exceptional cases and not for permanent use)

Protection against highspeed particles. Symbols for lens and frame:

F(T)	Low Energy Impact	45 m
B(T)	Medium Energy Impact	120 m
A(T)	High Energy Impact	190 m

If the F, B and A symbols are not the same for the visor and the headgear, then the lowest level should be assigned to the complete visor and face protection. The letter T immediately after the impact letter allows a use for high-speed particles at extremes of temperature. In the absence of the letter T, the protector must be used in ambient temperature.

CLEANING AND DISINFECTION: Face protector must be cleaned or disinfected only with a solution very low in alcohol on the headgear. For the visor, apply soft brush or air spray to remove any particles before transport or storage.**MAINTENANCE:** Like all other protective items, the life span depends on the type of application, care and maintenance. If the visor or headgear is damaged, replace the entire article. We recommend that all models are kept in dry and clean condition when not in use. In every case, the product must be replaced after up to a period of three years **REPLACEMENT VISOR:** Make sure that the spare visor are identical to the model no. of origin. **TRANSPORT:** The protector must be transported in the original packing. **STORAGE:** Store between 5 to 55°C; humidity less than 90%**Declaration of Conformity:** <https://www.hc-safety.com/files/EU-DoC-F03-F03A-2754-0538-085-12-20-1048-.pdf>**ESPAÑOL -MANUAL USO DE LA PANTALLA FACIAL**

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