
METRIA

Digital Chronometer, RS-021

Please read the User Manual carefully before use, and follow all operating and safety instructions!

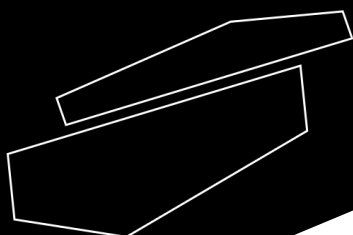user manual

english

User Manual

RS-021 Digital Chronometer

Stopwatch Function

In the normal running time state, press the C or D key (the C and D keys have the same function) to switch to stopwatch mode. If the stopwatch display is not at zero, press B to stop the stopwatch, and then press A to reset it to zero.

1. Stopwatch Timing


Press B to start timing, then press B again to stop timing. (Press B repeatedly to continue starting/stopping timing.) Press A to reset the stopwatch to zero.

2. Split Timing

Press B to start timing. Press A to display the first period's time, then press A twice quickly to display the second period's time. Press A twice quickly again to display the latest time. Press B to stop the timing. Press A to reset the stopwatch to zero.

3. Second Stage Timing

Press B to start timing. Press A to display the first period's time, then press B to stop timing. Press A to display the second period's time, and press A again to reset the stopwatch to zero.

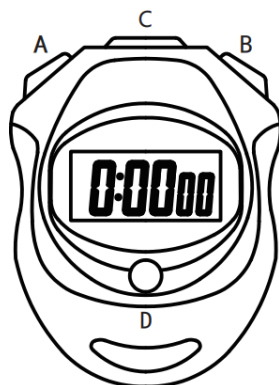Countdown Function

1. In time display mode, press C or D twice to enter countdown mode. The letters "TR" will appear in the upper left corner.
2. Press and hold A to enter the countdown setting (hour-minute-second sequence). Press B to adjust the time, or press and hold B to adjust quickly. Press C or D to confirm and complete the setting.
3. After setting, press B to start the countdown. Press B again to pause, and press it again to continue. Repeat this until the countdown finishes. When the countdown ends, a beep will sound for 30 seconds.
4. To reset after setting, press B to start the countdown, press B again to pause, and press A to return to the set countdown time. Press and hold A to reset the countdown to zero.

METRIA

Counter Function

1. In time display mode, press C or D three times to enter the counter mode.
2. Press B to start counting. Each time you press B, the count will increase by one.
3. Press and hold A to reset the counter to zero.

Alarm Setting Function

1. In time display mode, press C or D four times to enter the alarm setting mode.
2. Set the alarm hours and minutes sequentially. Press A to switch between hours and minutes, and press B to adjust the time. Press and hold B to adjust quickly.
3. Press C or D to confirm the alarm time.

Clock Setting Function

1. In time display mode, press C or D five times to enter the clock setting mode. The first setting is for 12/24-hour mode. Press B to select between 12-hour or 24-hour mode, then press A to confirm and enter the time setting.
2. To set the hour and minute, press A to switch between them. Press B to adjust the value, or press and hold B to adjust quickly. Press A to confirm and proceed to the year setting.
3. In the year setting, press A to switch between digits. Press B to adjust, or press and hold B to adjust quickly. Press A to confirm and proceed to the month-day setting.
4. For the month and day setting, press A to switch between month and day. Press B to adjust, or press and hold B to adjust quickly. Press C or D to complete the adjustment.

Alarm and Hourly Time Chime Switches

1. In time display mode, press and hold A, then press B. The alarm icon will appear to indicate that the alarm is on, and it will disappear when the alarm is off.
2. In time display mode, press and hold A, then press C. If the alarm icon is displayed, the hourly time chime is on. If the icon disappears, the hourly time chime is off.

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Important note for electronic devices sold in Spain

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Important note for electronic devices sold in France

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