



USER MANUAL

The 24h Clock (timer), reference TIME-W24-001, has a Clock function (12/24 hours), Stopwatch function and Countdown timer function.

NOTE: One LR44H button battery is included. Remove the protective paper from inside the battery compartment before using it for the first time.

I. 12/24 hour setting

Select the 12/24 system using the red selector on the side.

II. Time setting

- a. Press the MODE button to enter CLOCK mode.
- b. Adjust the hours, minutes and seconds by means of the buttons H, M, S respectively. By pressing each key for more than 2 seconds, the advance is faster.

III. Timer Operation

- a. Press the MODE button to enter TIMER mode, 00:00:00 appears on the display.
- b. Press the START button to start timing (COUNT-UP mode)
- c. To stop it, press the STOP button.
- d. Press the RESET button to reset the values.

IV. Countdown timer operation.

- a. Press the MODE button to enter TIMER mode, 00:00:00 appears on the display.
- b. Set the desired countdown time using buttons H, M, S respectively.
- c. Press the START button to start the countdown.
- d. If you want to pause the countdown, press the START button again.
- e. To stop it, press the STOP button.
- F. Press the RESET button to reset the values.

V. Memory function

When the time is up, the last entered time record can be accessed by pressing the MEMORY button.