
## R20 Chemical Resistant Vacuum Pump

*Please read the User Manual carefully before use, and follow all operating and safety instructions!*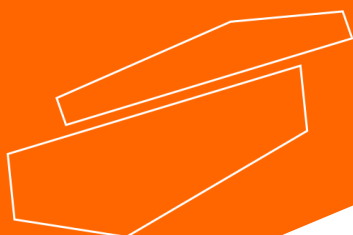

# user manual

english / español / français

# User Manual

## R20 series Chemically-resistant vacuum pump, 18 l/min, -750 mmHg

### Important notice

The vacuum pump is designed for laboratory use only. Before the setup and operation of the unit, please read these instructions carefully to familiarize you with all installation and operation process. Any individual should be well trained by reading the instructions or by technical people from the authorized distributors before operating the unit.

Do not modify or alter the unit in any way. Any modification or alteration will void the warranty and would result in potential hazards. We won't be responsible for any injury or damage caused by using the unit in case of any non-intended purpose, or modifying the unit by any person who is not authorized.

1. **When finish pumping, please do not turn off the pump at once and continue to run the vacuum pump for at least two minutes in order to draw out the mist and tiny liquid to prolong the service life of the pump.**
2. The filter cartridge (optional) is used to absorb moisture and dust. Please replace it when it is out of function to maintain a high pumping efficiency.
3. If the pump does not work when turning on the switch, please release the vacuum and turn-on again. If still fail to start the pump, there may be a problem in the switch or motor, please contact local distributor for help immediately.
4. Never use the pump with any flammable gas or toxic material.
5. In case of any damage with the pump, please directly contact local distributor or us for service. Do not repair it by yourself to avoid hazard.
6. If the supply cord is damaged, it must be replaced by the manufacturer, its service agent or similarly qualified persons in order to avoid a hazard.

## Service

In order to guarantee this equipment Works safely and efficiently, it must have a regular maintenance. In case of any faults, do not try to repair it yourself. If help is needed, you can always contact your dealer or Labbox via [www.labbox.com](http://www.labbox.com)


Please provide the customer care representative with the following information:

- Serial number
- Description of problem
- Your contact information

## Installation

1. Check voltage specified on name-plate on the bottom. Make sure if it matches the line voltage in your location.
2. Install the instrument in a clean, dust-less and ventilated area under 40°C.
3. To enable suction function, connect the inlet of the pump to the outlet of objective equipment with a high-pressure hose.

## Assembly & Disassembly diagram

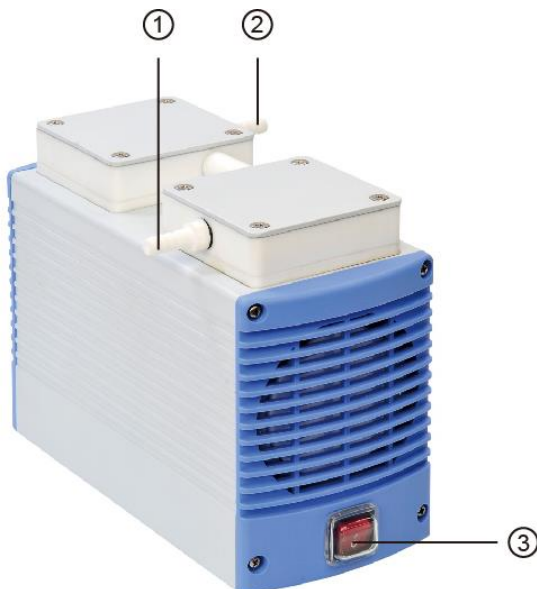Position	Designation
1	Air inlet
2	Air outlet
3	On/Off switch

# Operation

## Operating procedure

1. The diagram of the pump system is shown as above.
2. Once all tubes have been connected to Inlet (1) Outlet (2) properly then activate the pumping function by pressing (3) on/off button.
3. To adjust vacuum of the pump, an optional vacuum regulator/ moisture trap is needed.

**Warning:** When pump is used in water filtration, please do not make the liquid level of flask exceed safety level. Failure to comply could result in serious damage to the pump and void the warranty. If it occurs, please contact the distributor immediately for service.

## Setup moisture trap (Option)

Moisture trap (Vacuum regulator) is an optional item that is used to adjust vacuum and absorb moisture and particle.

The flow chart of installing vacuum regulator/moisture trap

	Remove the Teflon tier connector at side of cylinder head with a spanner.
	Fix the moisture trap at the location of the Teflon tier connector removed from Step 1.
	Fix the glass bottle. Finished

## Warranty

This instrument is warranted to be free from defects in materials and workmanship under normal use and service, for a period of 48 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claim under the warranty please contact your supplier.

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